




Cadenza CD-C18

75 x 3 mm

Application

**Volicitin-related Compounds  
from the Oral Secretion of Caterpillars**

鱗翅目昆虫幼虫の唾液中ポリシチン関連化合物

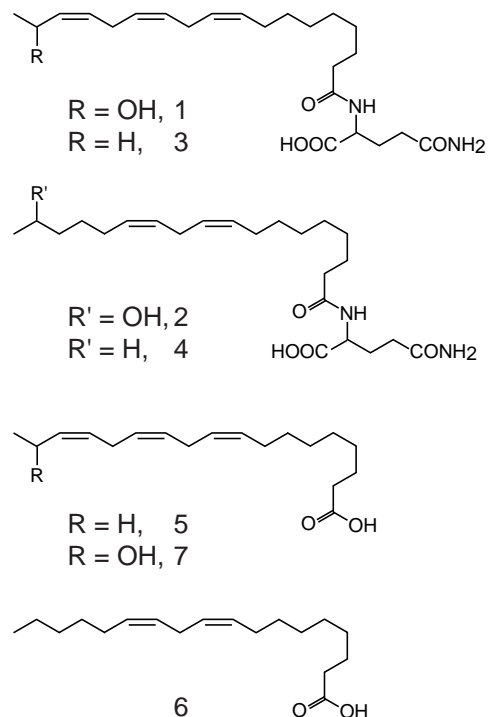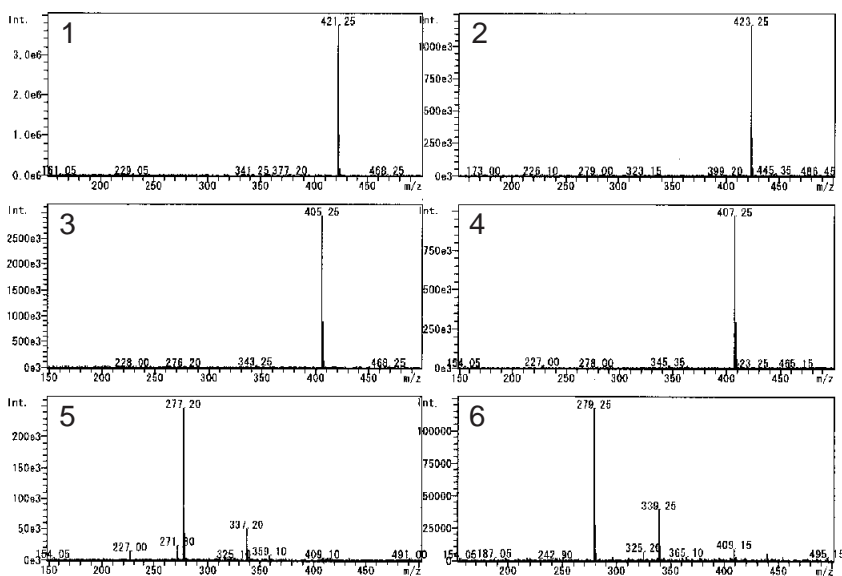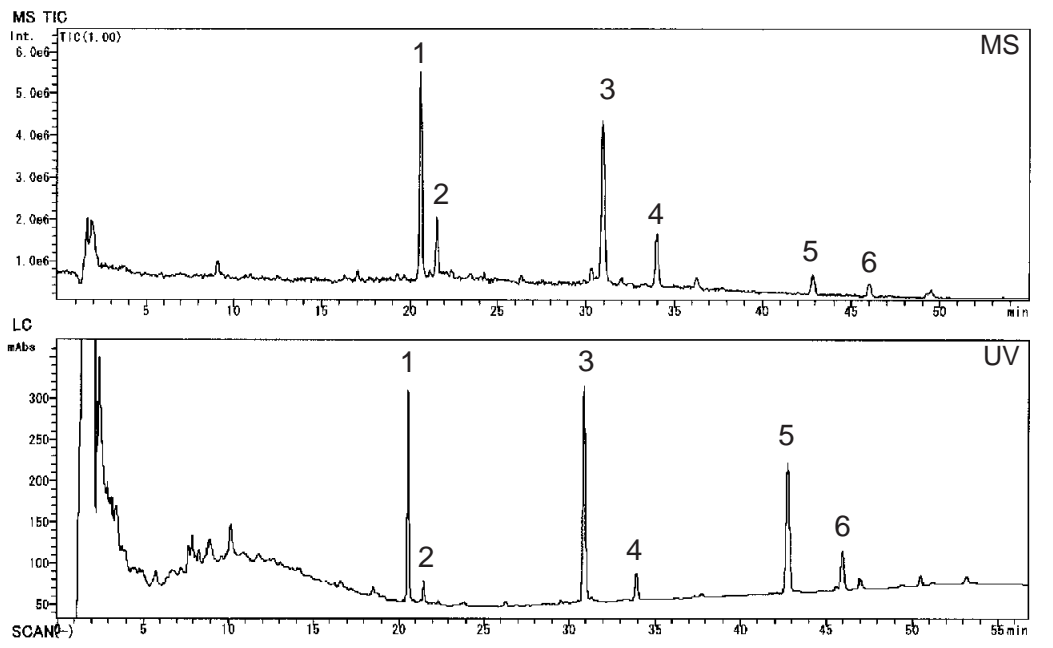

Cadenza CD-C18, 75 x 3 mm  
 A: 0.05% acetic acid in water  
 B: 0.08% acetic acid in acetonitrile  
 20-95%B (0-40min)  
 0.2 mL/min, 60°C, 10 μL  
 Negative ESI mass, UV at 200 nm

Courtesy of Dr. Naoki Mori, Kyoto Univ.