

Cadenza CD-C18


150 x 2 mm

Application


LC-MS/MS analysis of clenbuterol (CBR)

クレンブテロール (CBR) のLC-MS/MS分析(告示試験法)

CBR CBRd9 0.5ppb 0.5ppb STD imtakt

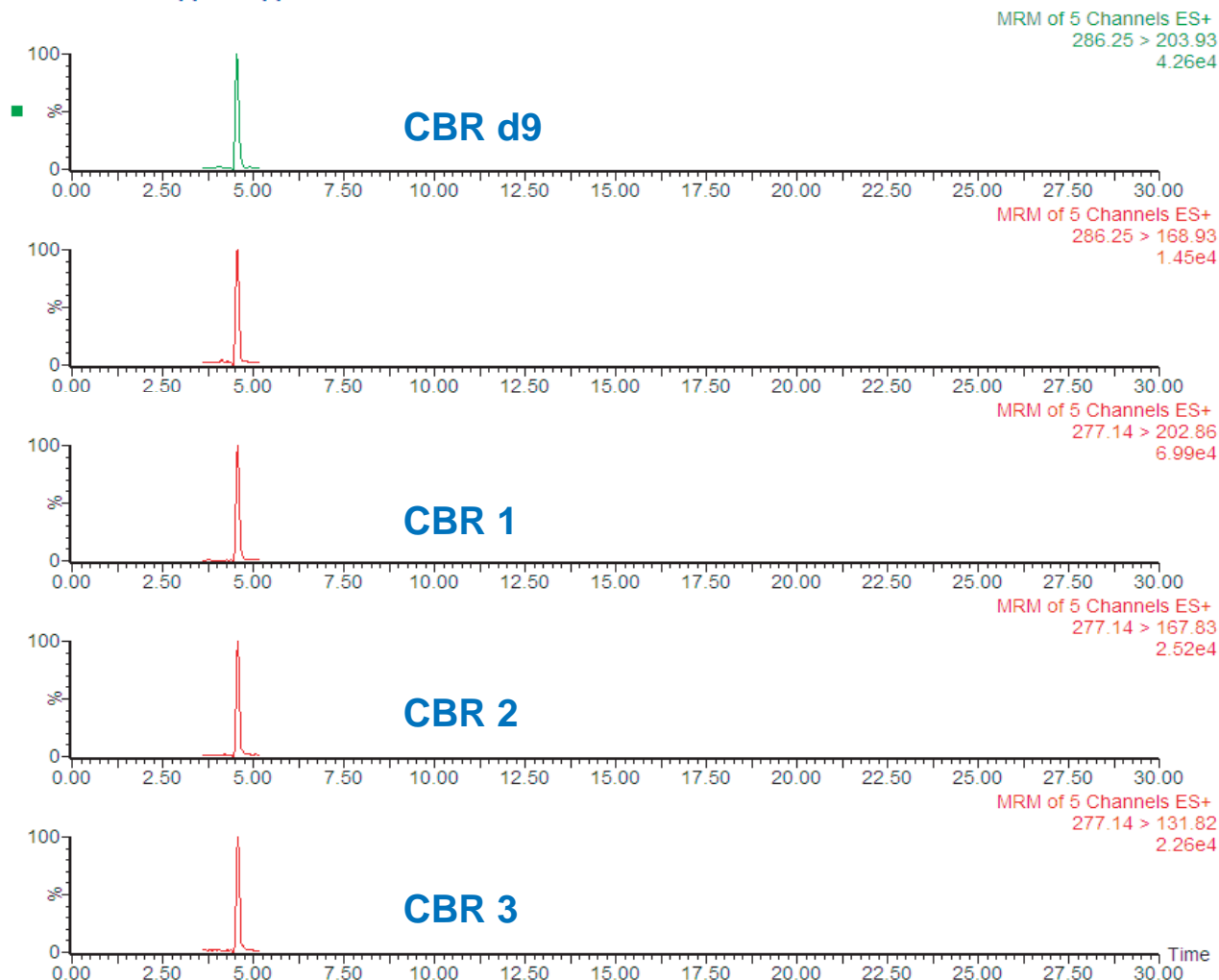

Cadenza CD-C18, 150 x 2 mm
 A: 0.1% formic acid in water
 B: 0.1% formic acid in acetonitrile
 5-95%B (0-7min), 95%B (7-10min)
 Runtime 15min
 0.2 mL/min, 40°C
 5uL (CBR 0.5ppb, IS 0.5ppb)
 ESI, MRM, positive

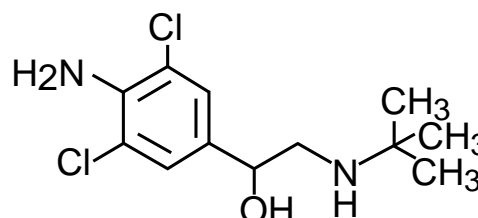

clenbuterol (CBR)

Courtesy of MAEDA Naoyuki, ASSOCIATION OF MEAT SCIENCE & TECHNOLOGY INSTITUTE, JAPAN