


Scherzo SM-C18
Unison UK-C18


75 x 3 mm

Application


Spermine and spermidine
スperlミンとスperlミジン

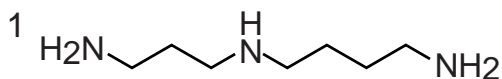

spermidine

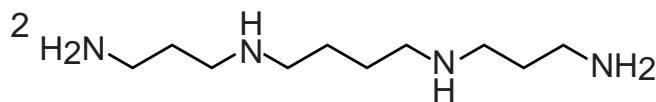

spermine

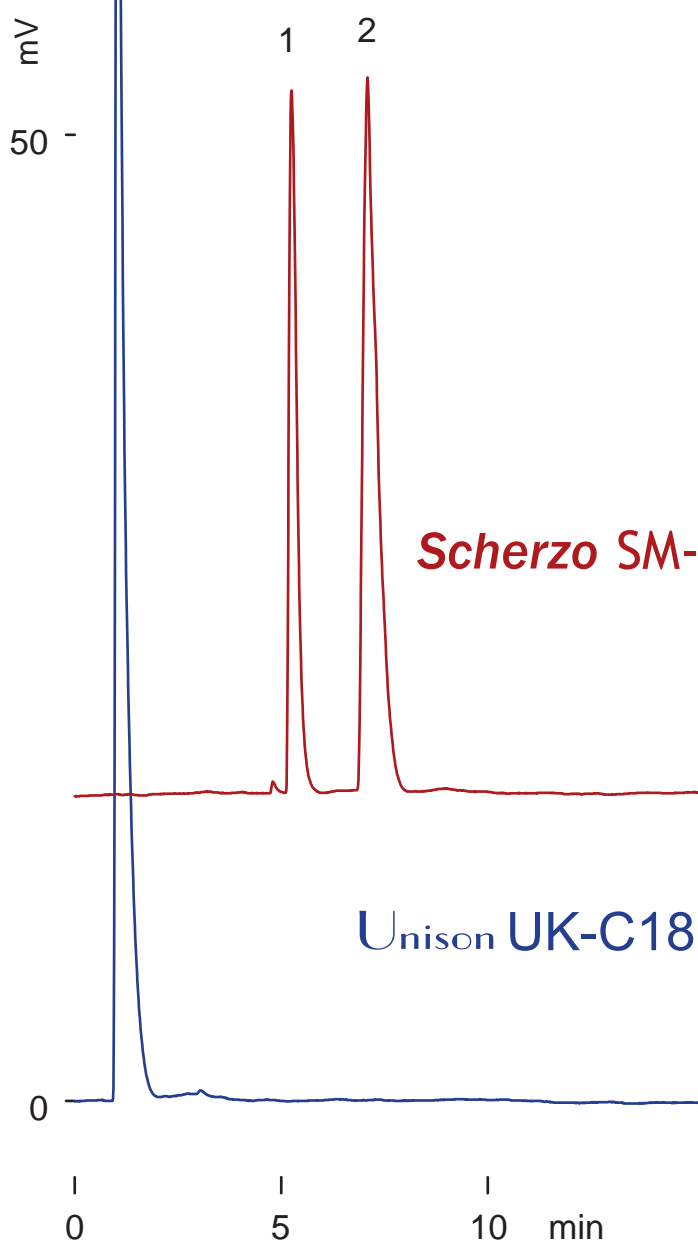

Scherzo SM-C18

Unison UK-C18

75 x 3 mm
 A: 10mM ammonium acetate
 B: 100mM acetic acid /acetonitrile
 = 50 /50
 0-40 %B (0-10min)
 0.4 mL/min (4MPa), 37 deg.C
 ELSD (spray chamber 10 deg.C,
 drift tube 45 deg.C)
 2 uL (1-2 ug)