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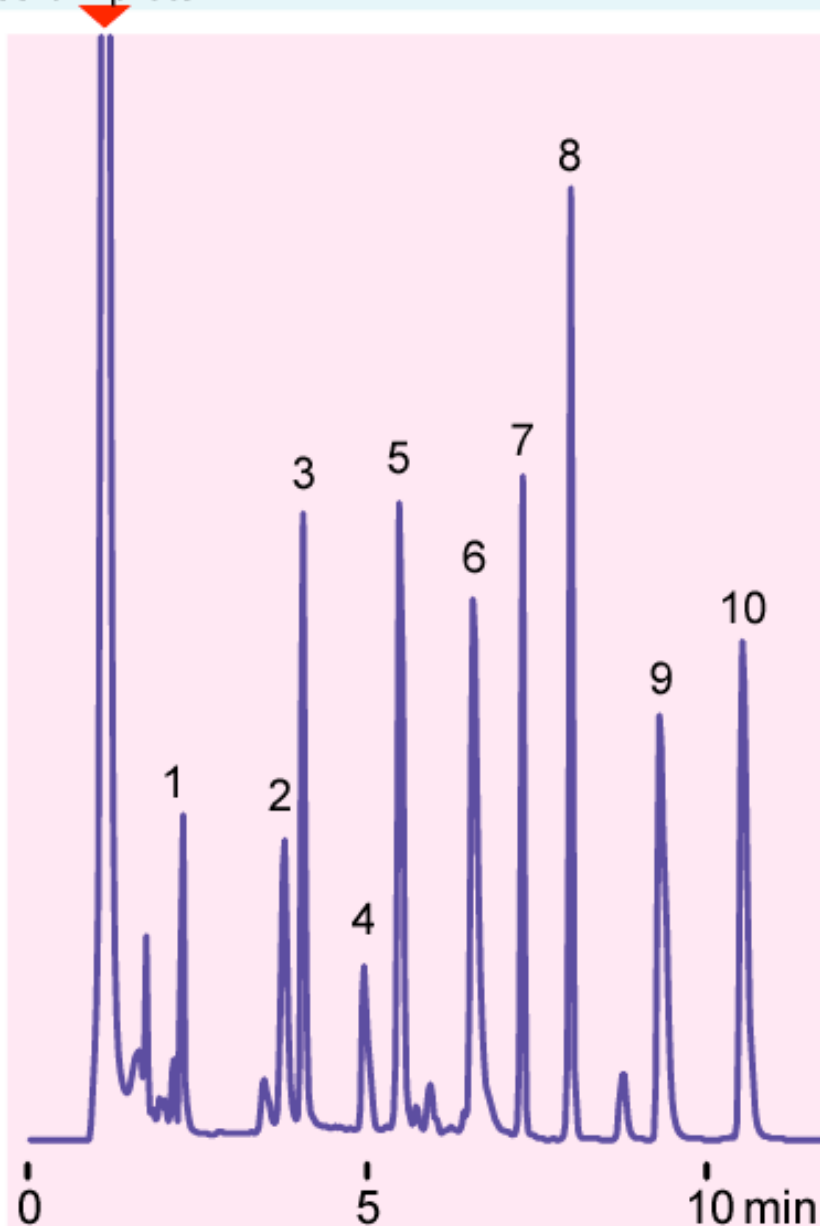
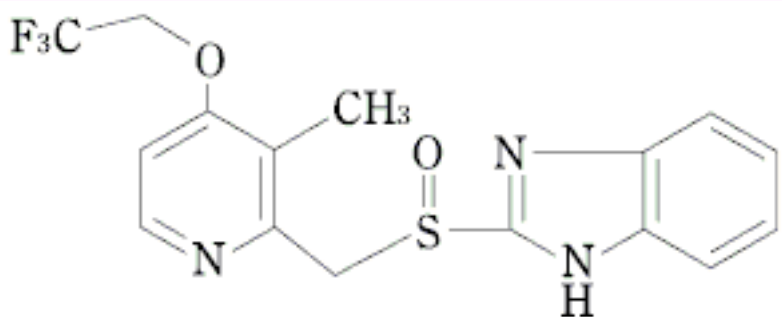
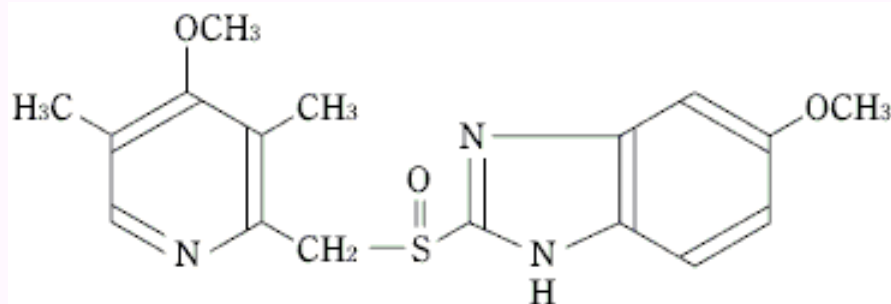
Omeprazole
(Internal Standard)

Sample Preparation

*Direct Injection vs.
Protein Precipitation*

Direct Injection

↓ Plasma
↓ Centrifugation
↓ Addition of I.S.
 9:1 H₂O (I.S.) to Plasma
↓ Injection

Protein Precipitation

↓ Plasma
↓ Centrifugation
↓ Protein Precipitation
 4:1 ACN (IS) to Plasma
 ↓ Filtration
↓ Injection

Analytical Conditions

*Direct Injection vs.
Protein Precipitation*

Direct Injection

Column

Cadenza HS-C18
20×4.6 mm, 3 µm (Imtakt)
with Guard Column

Mobile Phase

A, 10 mM ammonium formate
(pH3.0)

B, Acetonitrile

100 % A (0→0.5 min.)

0→100 % B (0.5→2.5 min.)

100 %B (2.5→3.5 min.)

Flow Rate 500 µL/min.

Temperature 40 deg. C

Injection Volume 10 µL

MS/MS ESI, Positive, MRM

Protein Precipitation

Column

ODS Column
20×4.6 mm, 5 µm
without Guard Column

Isocratic

A, 10 mM ammonium formate
(pH3.0)

B, Methanol

C, Acetonitrile

A/B/C = (27.5 : 35.0 : 37.5)

0 - 8.5 min.

Flow Rate 70 µL/min.

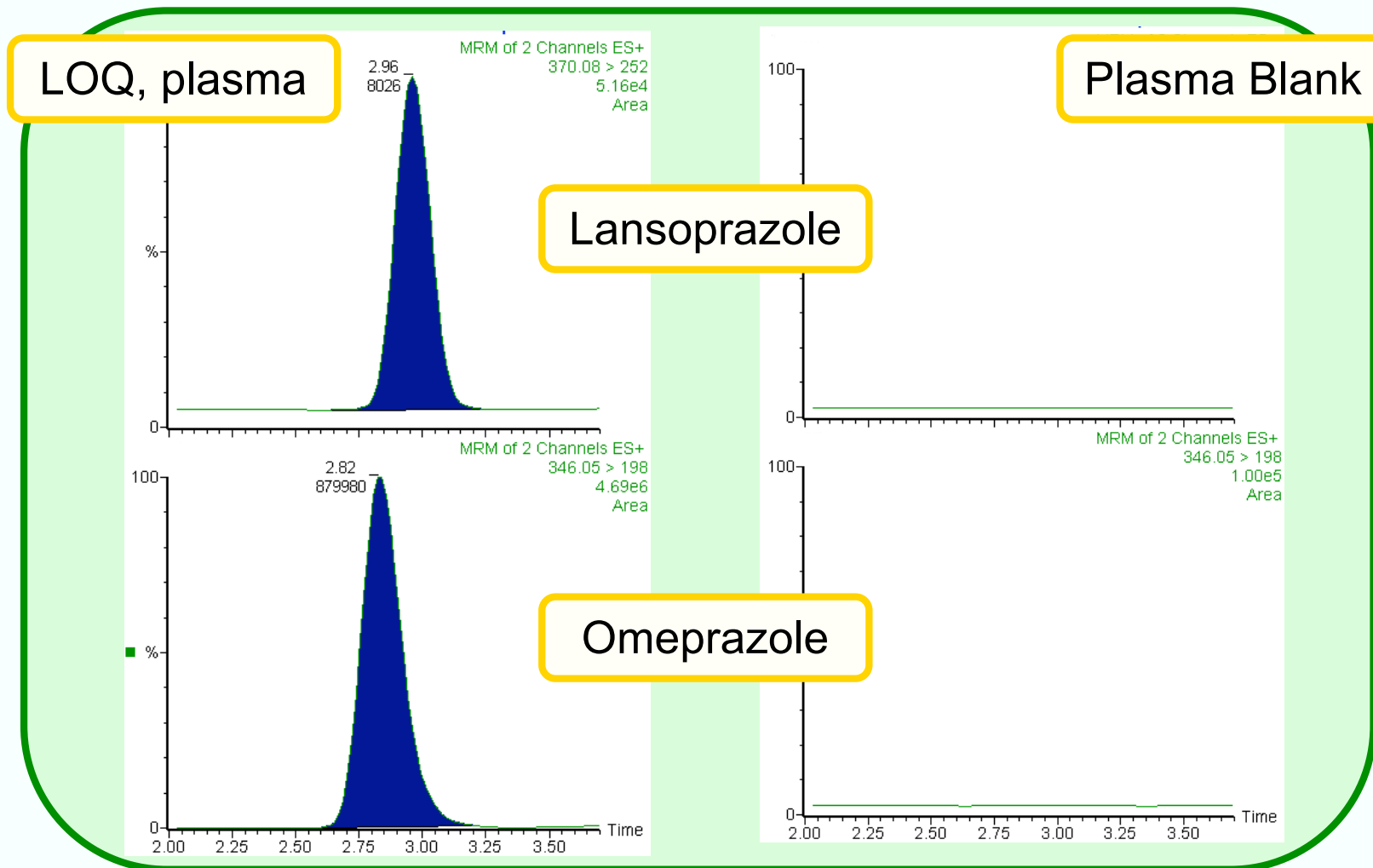
Temperature 40 deg. C

Injection Volume 1 µL

MS/MS ESI, Positive, MRM

Chromatogram


*Direct Injection vs.
Protein Precipitation*

TOWA PHARMACEUTICAL CO., LTD.

Linearity

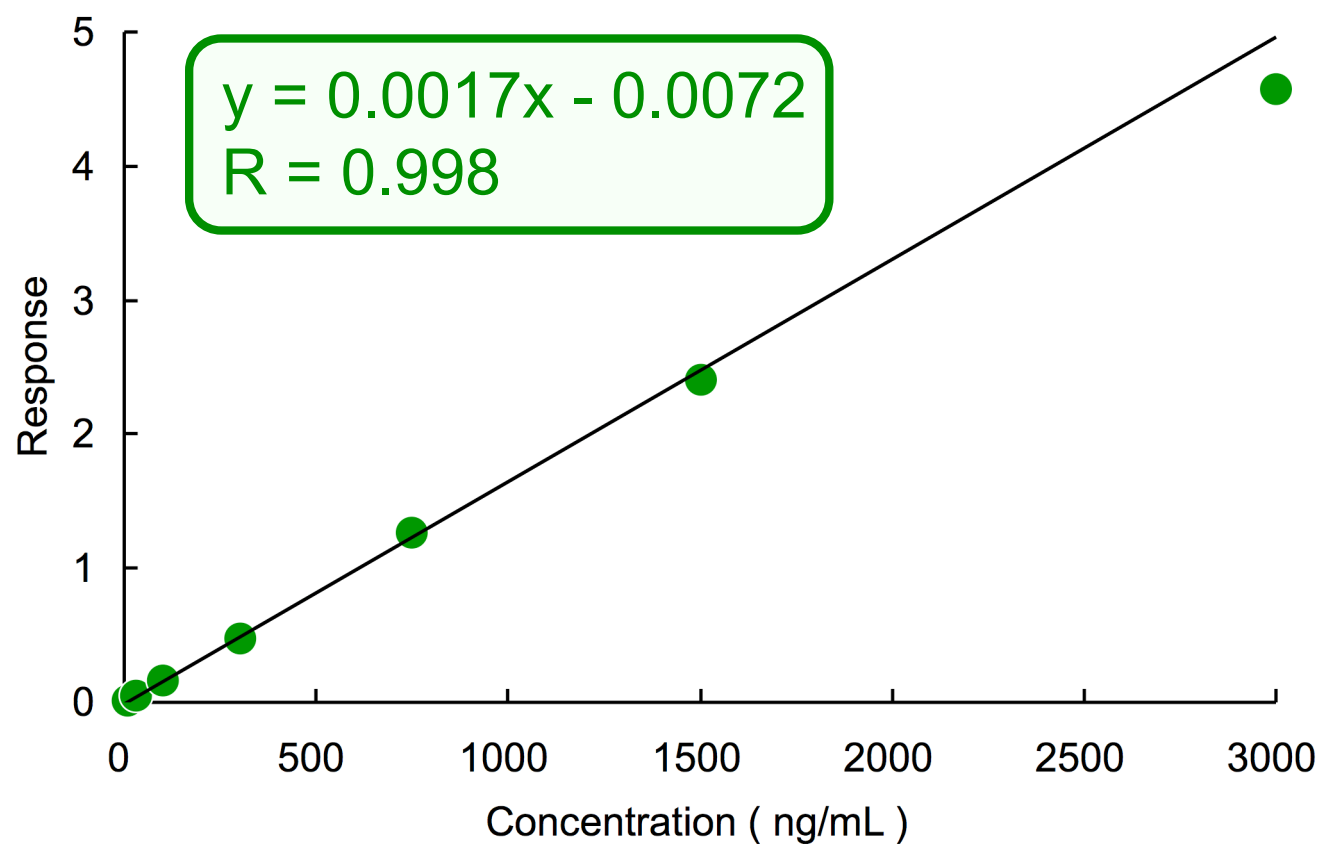*Direct Injection vs.
Protein Precipitation*



10 - 3000 ng/mL, Weight $1/x^2$, $n = 3$

TOWA PHARMACEUTICAL CO., LTD.

Linearity

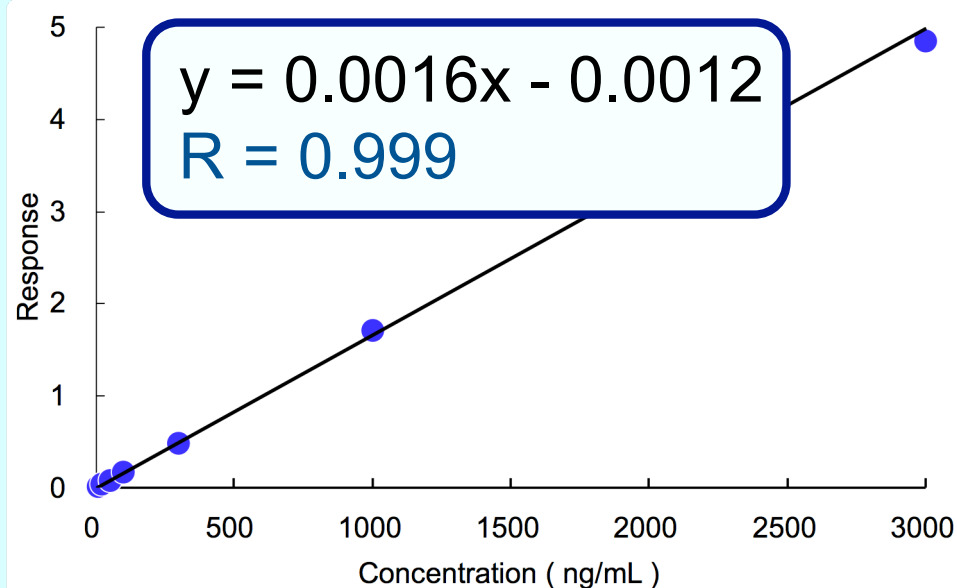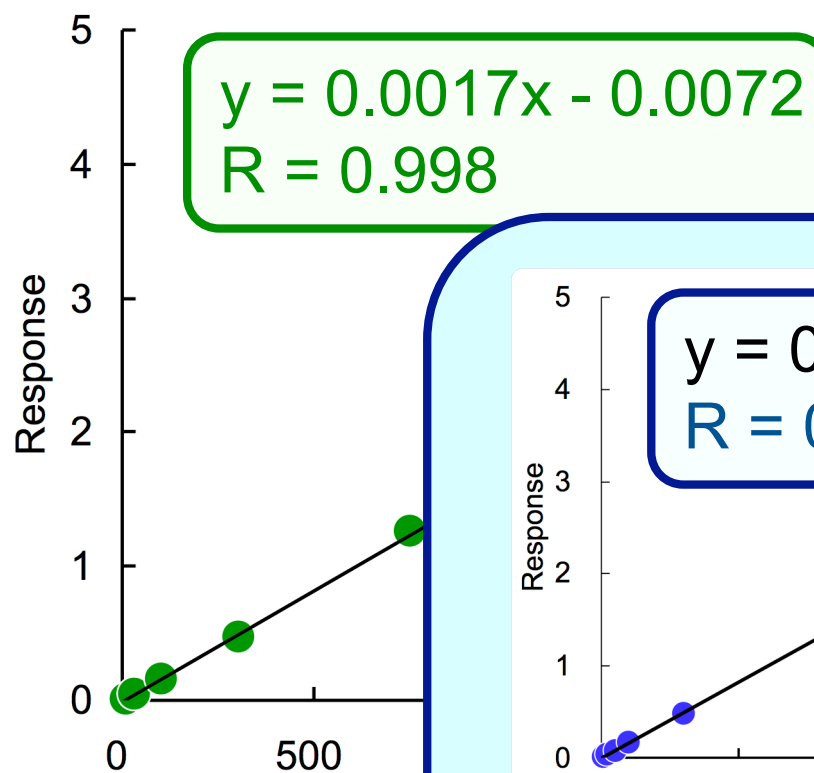*Direct Injection vs.
Protein Precipitation*

10 - 3000 ng/mL, Weight $1/x^2$, $n = 3$

TOWA PHARMACEUTICAL CO., LTD.

Accuracy and Precision


*Direct Injection vs.
Protein Precipitation*

	Nominal Concentration (ng/mL)					
	10		300		3000	
Accuracy (Mean, %)	-1.9	5.5	4.3	-0.4	-6.4	0.7
Precision (CV, %)						
Intra-batch (n = 3)	2.7	4.0	2.4	3.1	2.5	2.8
Inter-batch (n = 6)	5.0	6.1	4.0	5.1	6.6	4.8

Quality Control

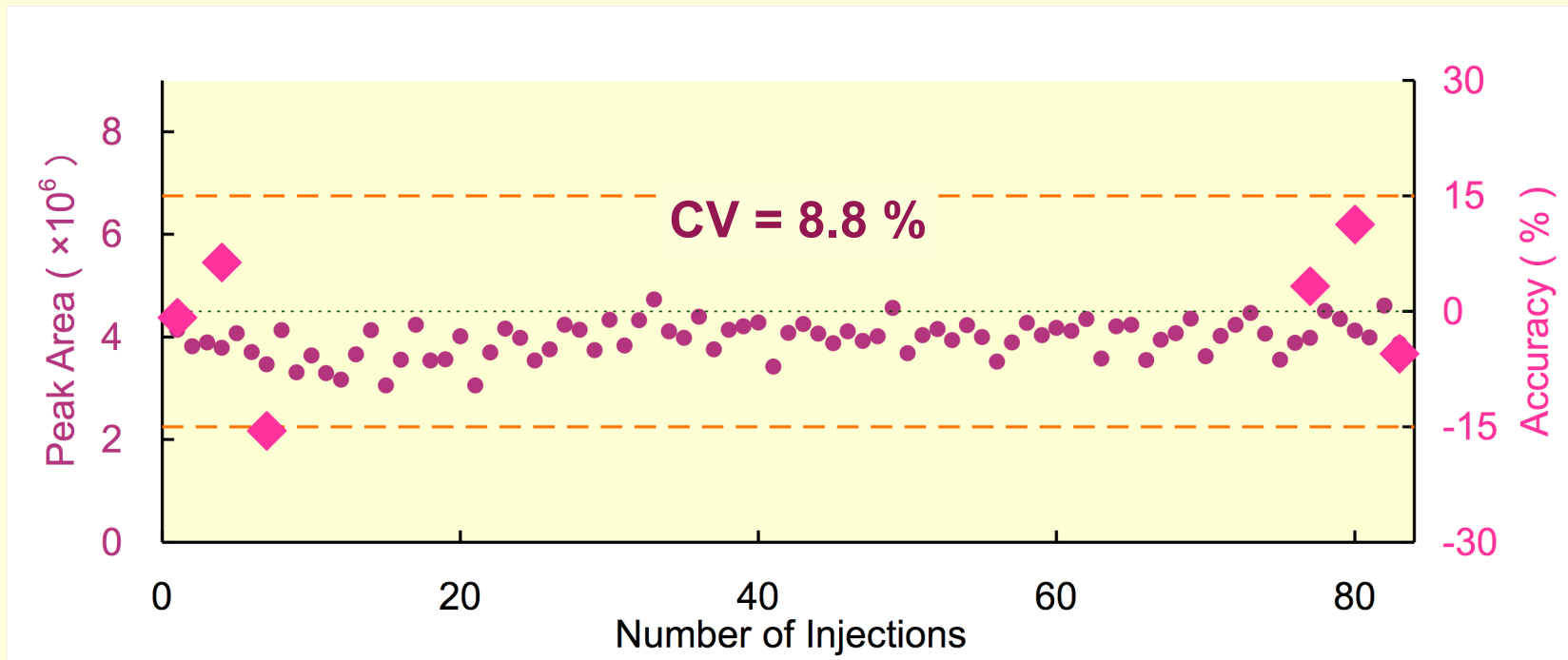Clinical Trial

Lansoprazole 30 mg Single Dosa (One Healthy Adult male)

● - IS Peak Area

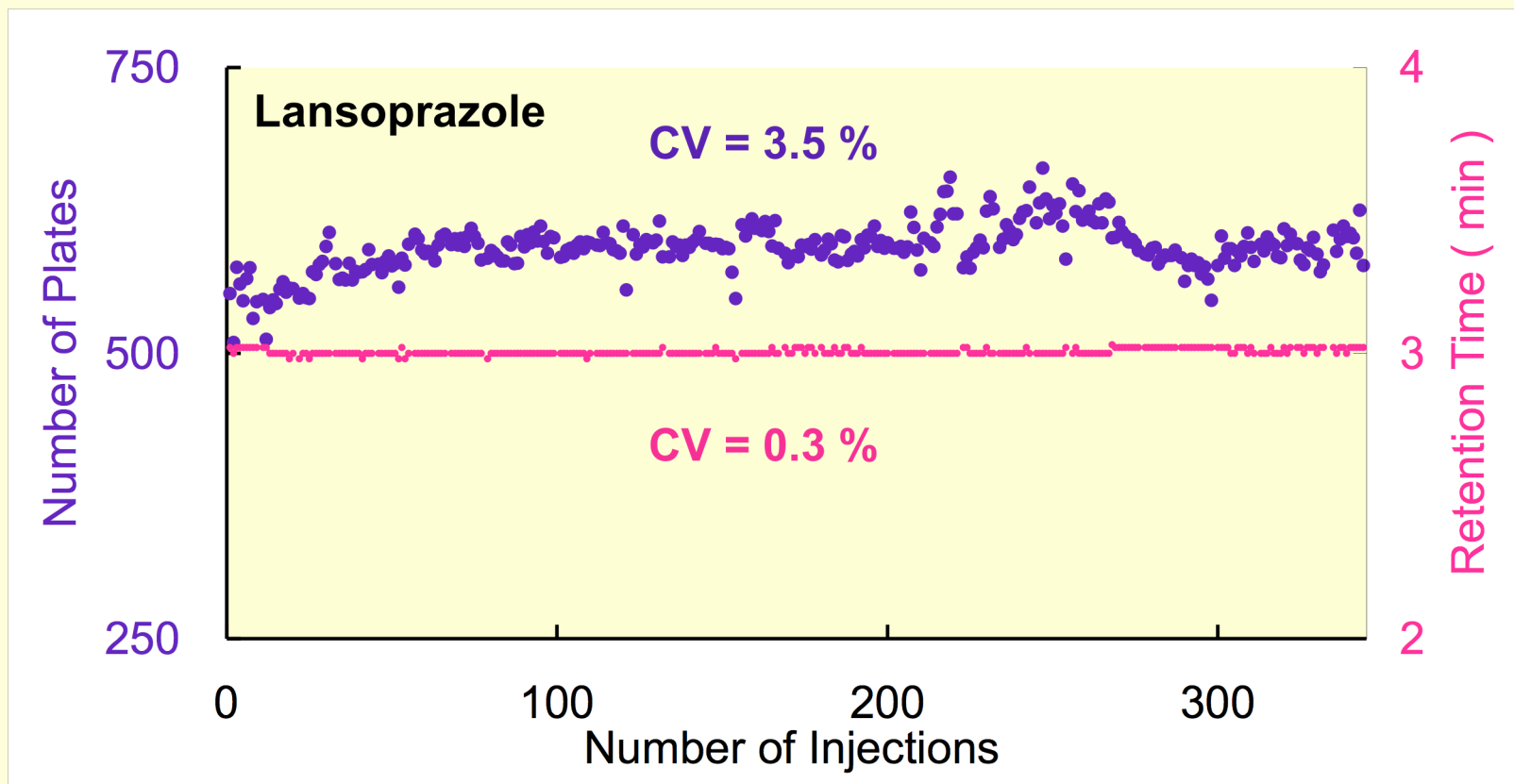
◆ - QC Samples (10, 300 and 3000 ng/mL); QC Range = ± 15 %

TOWA PHARMACEUTICAL CO., LTD.

Longevity


Clinical Trial

TOWA PHARMACEUTICAL CO., LTD.

Plasma Concentrations

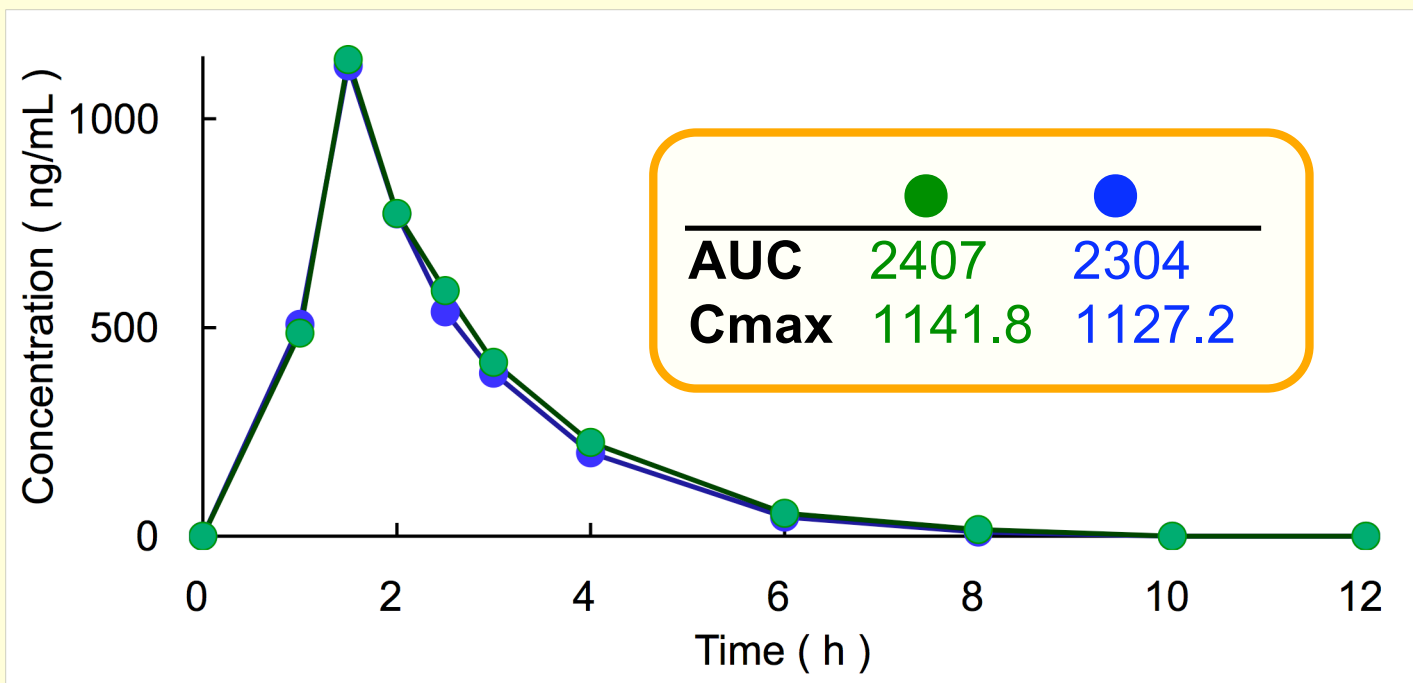Clinical Trial

- - Direct Injection
- - Protein Precipitation

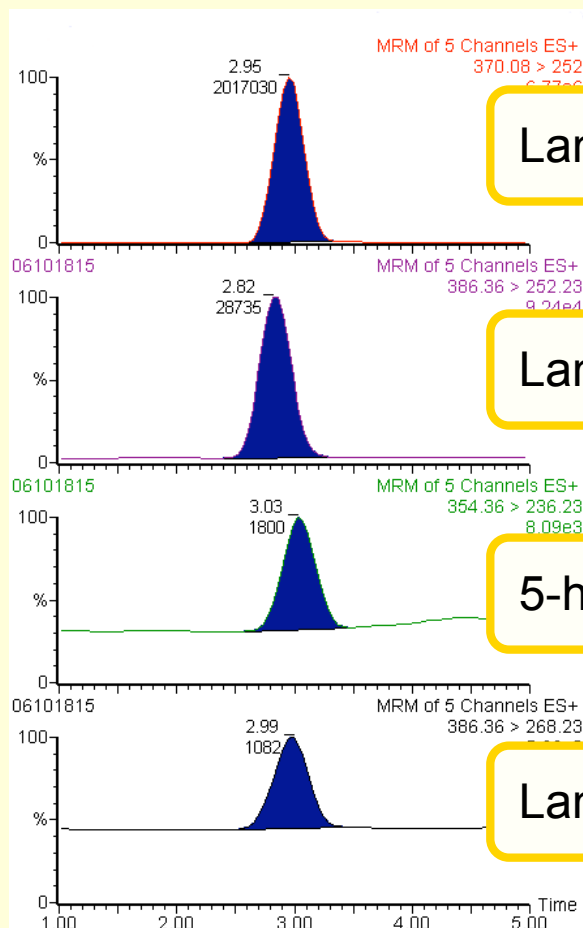
AUC - ng·h/mL
Cmax - ng/mL

TOWA PHARMACEUTICAL CO., LTD.

Metabolites

Clinical Trial

Lansoprazole

Lansoprazole Sulfide

5-hydroxy Lansoprazole

Lansoprazole Sulfone

TOWA PHARMACEUTICAL CO., LTD.

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**Novel Hydrophobic/Hydrophilic Hybrid ODS Column
Cadenza HS-C18**

**Direct Injection LC-MS/MS Analysis of
Lansoprazole in human plasma**

**Method Validation
Clinical Trial
Direct Injection vs. Protein precipitation**

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**Novel Hydrophobic/Hydrophilic Hybrid ODS Column
Cadenza HS-C18**

**Direct Injection LC-MS/MS Analysis of
Lansoprazole in human plasma**

Satisfactory

Direct injection vs. Protein precipitation

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**Novel Hydrophobic/Hydrophilic Hybrid ODS Column
Cadenza HS-C18**

**Direct Injection LC-MS/MS Analysis of
Lansoprazole in human plasma**

Direct Injection vs. Precipitation

Excellent?

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**Novel Hydrophobic/Hydrophilic Hybrid ODS Column
Cadenza HS-C18**

**Direct Injection LC-MS/MS Analysis of
Lansoprazole in human plasma**

Direct Injection vs. Extraction

TOWA PHARMACEUTICAL CO., LTD.