



Unison UK-Amino
Unison UK-Silica

150 x 2 mm

Application


PQQ: pyrroloquinoline quinone (Normal Phase)

新ビタミン PQQ:ピロロキノリンキノンの順相分析

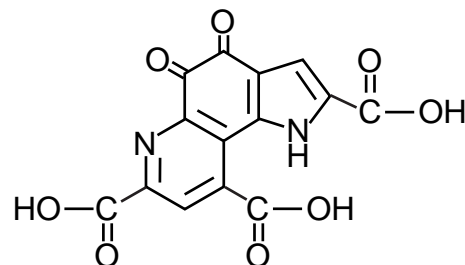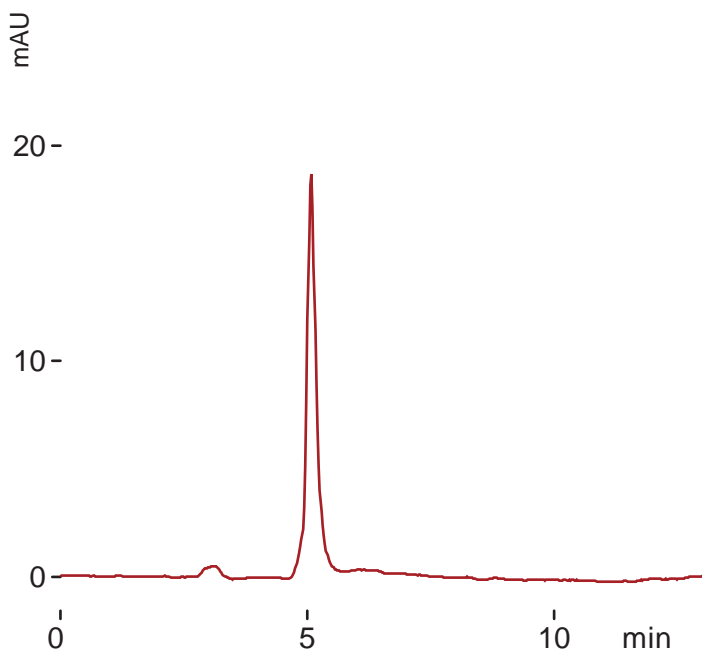

PQQ (pyrroloquinoline quinone)

Unison UK-Amino, 150 x 2 mm
A: acetone / formic acid = 100 / 0.1
B: water / formic acid = 100 / 0.1
0 - 100%B (0-10min)
0.2 mL/min (3MPa)
37 deg.C, ELSD, 2 uL (2ug)

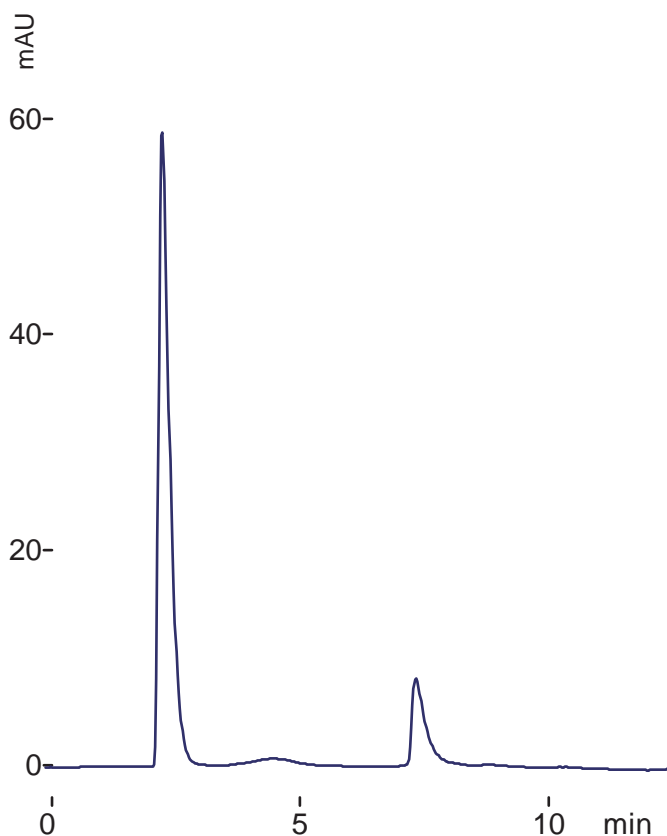

Unison UK-Silica, 150 x 2 mm
A: acetone / formic acid = 100 / 0.1
B: water / formic acid = 100 / 0.1
0 - 100%B (0-10min)
0.2 mL/min (3MPa)
37 deg.C, ELSD, 2 uL (2ug)