


Unison UK-Amino

250 x 3 mm


Application

Unsaturated Chondro-Disaccharides


不飽和コンドロ二糖

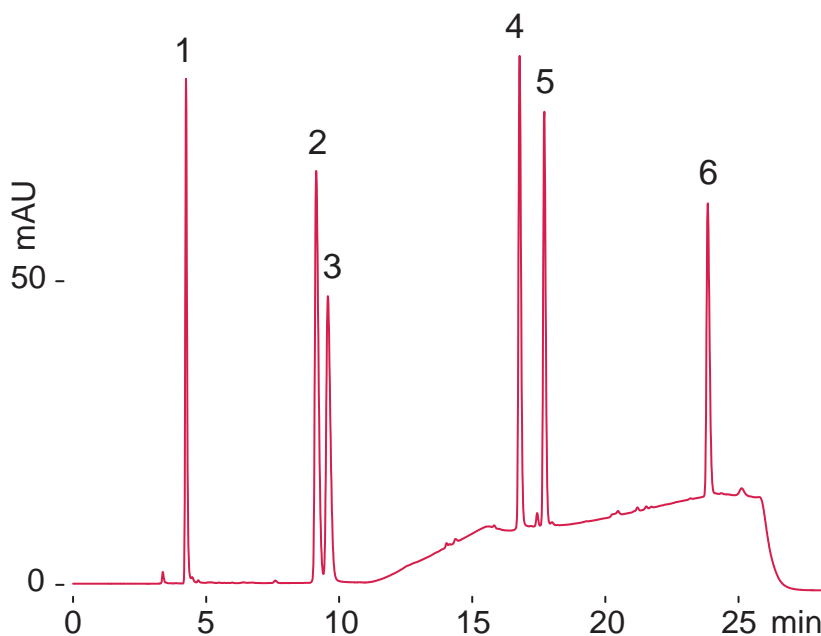


Unison UK-Amino, 250 x 3 mm
 A: 10mM NH₄HCO₃, B: 400mM NH₄HCO₃
 10%B(0-7min), 10-100%B(7-20min), 100%B(20-22min)
 0.4 mL/min(13MPa), 50 deg.C, 232 nm, 5 uL(0.4-1.5 ug)

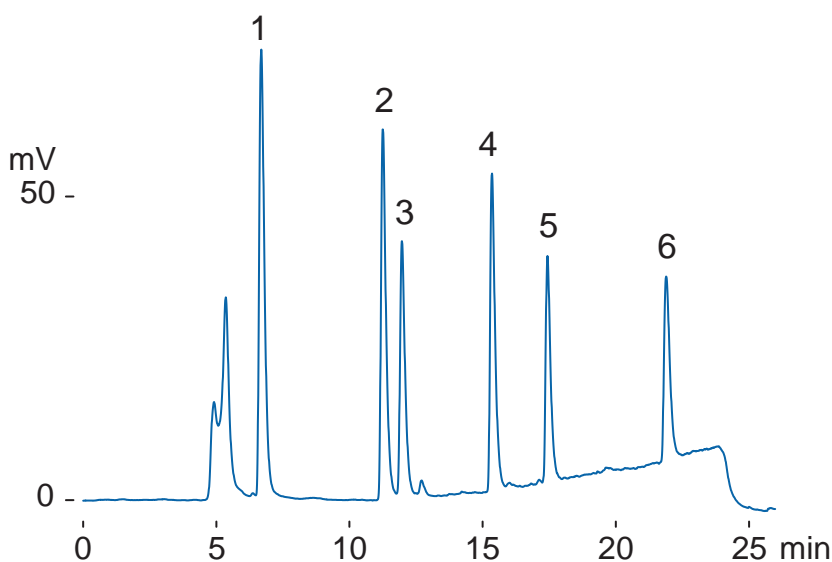


Unison UK-Amino, 250 x 3 mm
 A: methanol, B: water /TEA /AcOH = 1000 /41 /16
 3-90%B (0-20min)
 0.4 mL/min (8MPa), 50 deg.C, ELSD, 10 uL (1-3ug)

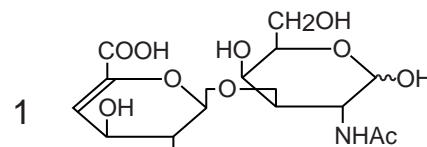


Δ Di-0S

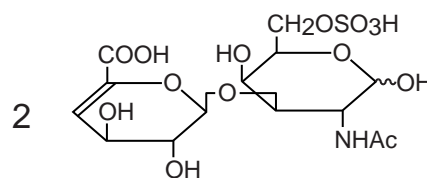


Δ Di-6S

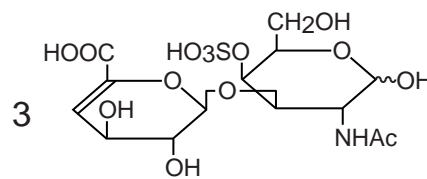


Δ Di-4S

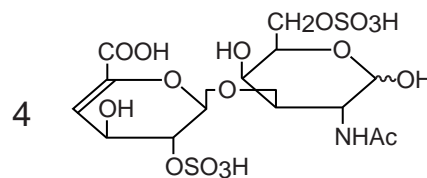

Δ Di-diS_D

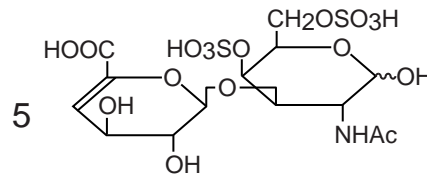

Δ Di-diS_E

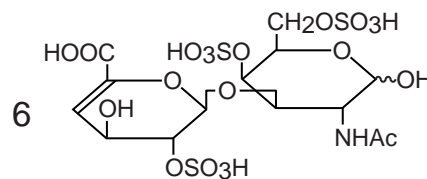

Δ Di-triS

Courtesy of Seikagaku Corporation, Japan