





Scherzo SM-C18
Unison UK-C18

30 x 3 mm

Application

Adenosine phosphates (AMP, ADP, ATP)

アデノシンリン酸 (AMP, ADP, ATP)

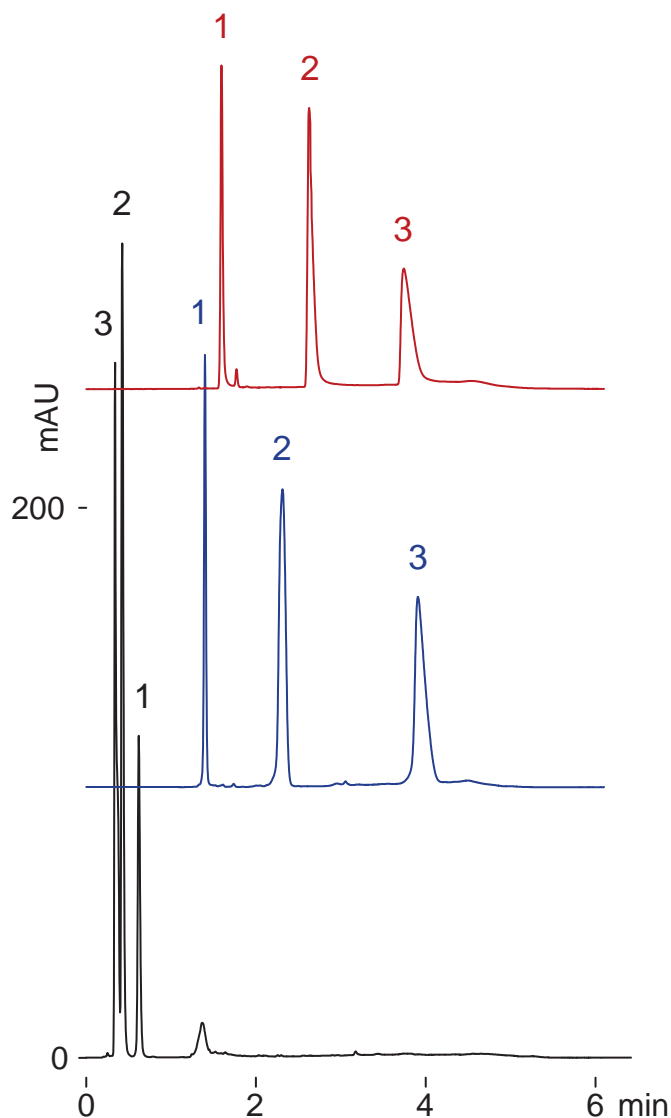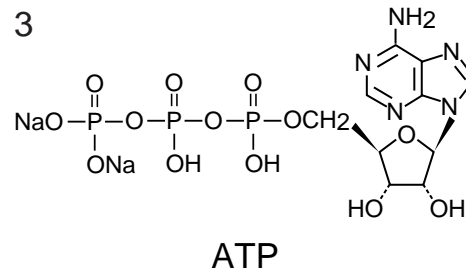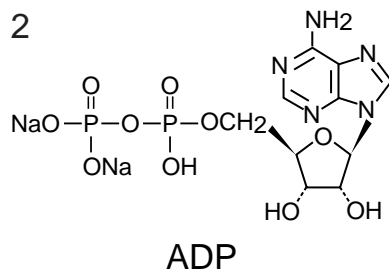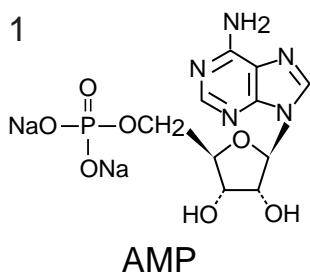

Scherzo SM-C18

A: 20mM CH₃COONH₄

B: 200mM CH₃COONH₄ /ACN = 70 /30

Scherzo SM-C18

A: 5mM NH₄H₂PO₄ : 5mM (NH₄)₂HPO₄
= 1:1, pH6.8

B: (20mM NH₄H₂PO₄ : 20mM (NH₄)₂HPO₄
= 1:1, pH6.8) /ACN = 50 /50

Unison UK-C18

A: 5mM NH₄H₂PO₄ : 5mM (NH₄)₂HPO₄
= 1:1, pH6.8

B: (20mM NH₄H₂PO₄ : 20mM (NH₄)₂HPO₄
= 1:1, pH6.8) /ACN = 50 /50

30 x 3 mm

0-100 %B (0-10min)

0.6 mL/min (5-6 MPa), 37deg.C, 260 nm, 0.8 uL (0.1-0.3ug)