


Cadenza CW-C18
Unison **UK-Phenyl**

75 x 3 mm


Application

Phthalate esters


フタル酸エステル類

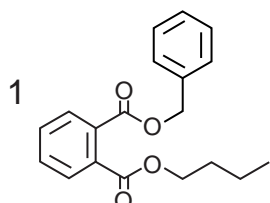


butylbenzyl phthalate
BBP

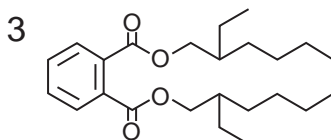


di-2-ethylhexyl phthalate
DEHP

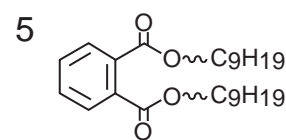


diisononyl phthalate
DINP

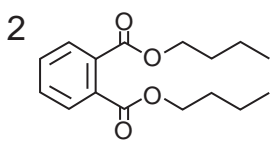


dibutyl phthalate
DBP

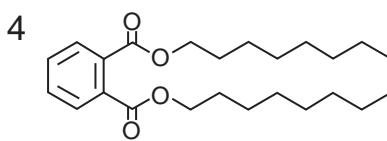

di-n-octyl phthalate
DNOP

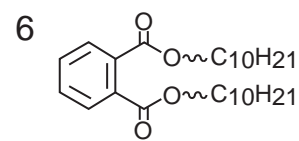

diisodecyl phthalate
DIDP

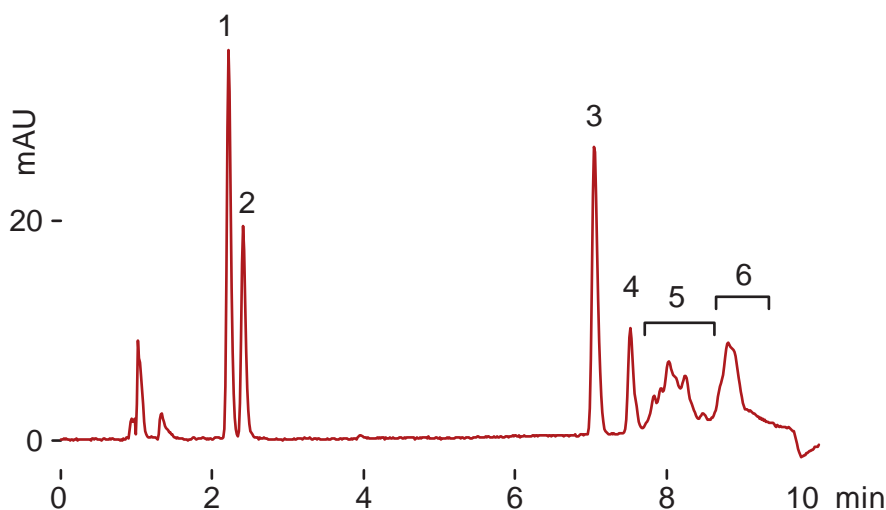

Cadenza CW-C18

75 x 3 mm

A: water / acetic acid
= 100 / 0.1

B: acetonitrile / acetic acid
= 100 / 0.1


70-100 %B (0-8min)

70 %B (8-10min)

0.4 mL/min (4MPa)

37 deg.C, 275 nm

2 uL (0.1-0.5ug)

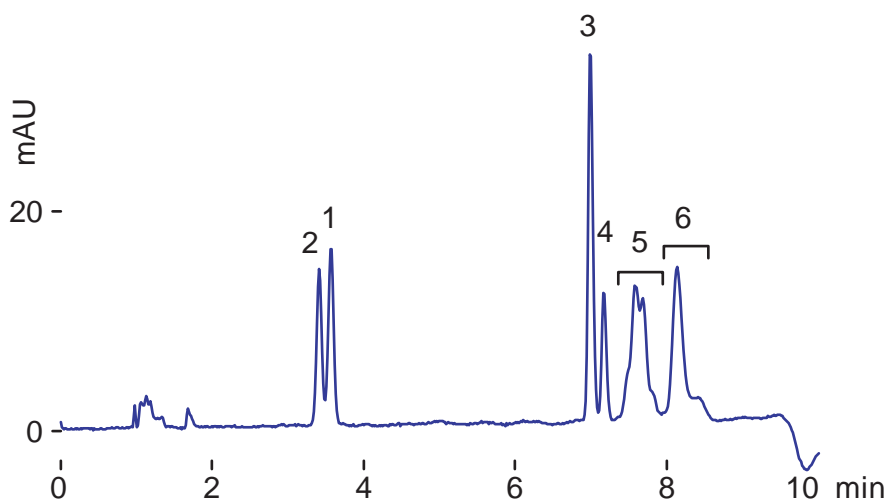

Unison UK-Phenyl

75 x 3 mm

A: water / acetic acid
= 100 / 0.1

B: methanol / acetic acid
= 100 / 0.1

70-90 %B (0-8min)

70 %B (8-10min)

0.4 mL/min (7MPa)

37 deg.C, 275 nm

2 uL (0.1-0.5ug)