


Presto FF-C18


50 x 4.6 mm
150 x 4.6 mm

Application

Pectin ペクチン

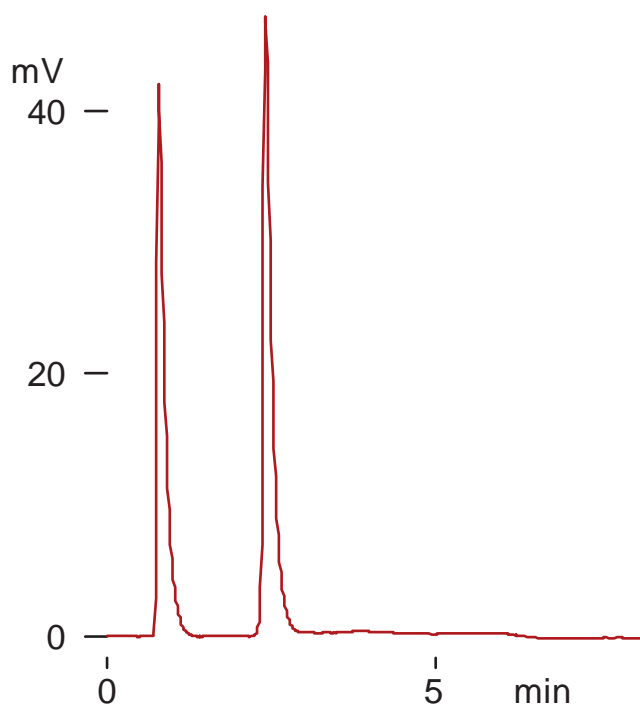

Presto FF-C18, 50 x 4.6 mm
A: water / formic acid = 100 / 0.1
B: acetonitrile / formic acid = 100 / 0.1
0-70 %B (0-5 min)
0.5 mL/min (7 MPa), 37 deg.C
ELSD (spray chamber 30 deg.C,
drift tube 80 deg.C)
2 uL (4 ug)

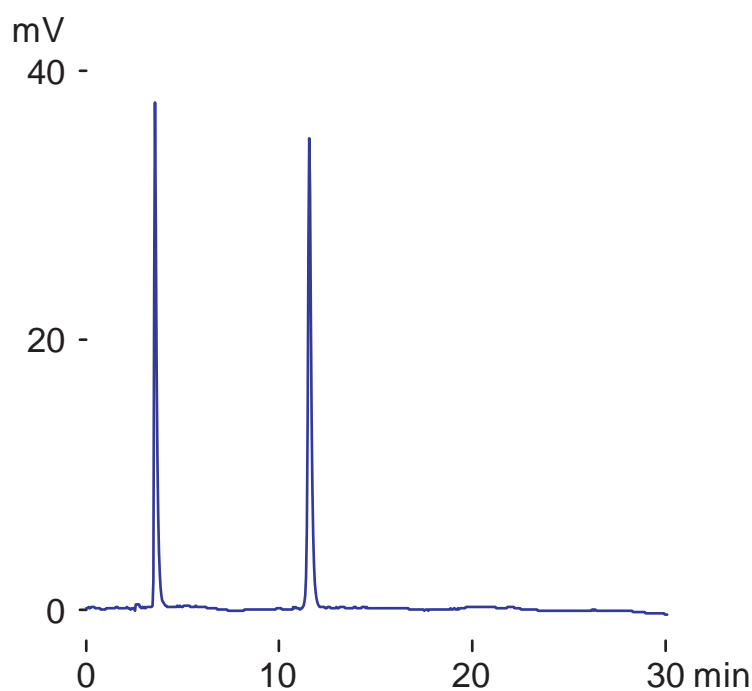

Presto FF-C18, 150 x 4.6 mm
A: water / formic acid = 100 / 0.1
B: acetonitrile / formic acid = 100 / 0.1
0-50 %B (0-25 min)
0.3 mL/min (11 MPa), 37 deg.C
ELSD (spray chamber 30 deg.C,
drift tube 80 deg.C)
2 uL (4 ug)