


Unison UK-Amino

250 x 3 mm

Application

High sensitivity analysis for saccharides using Corona CAD

Corona CAD による糖類の高感度分析

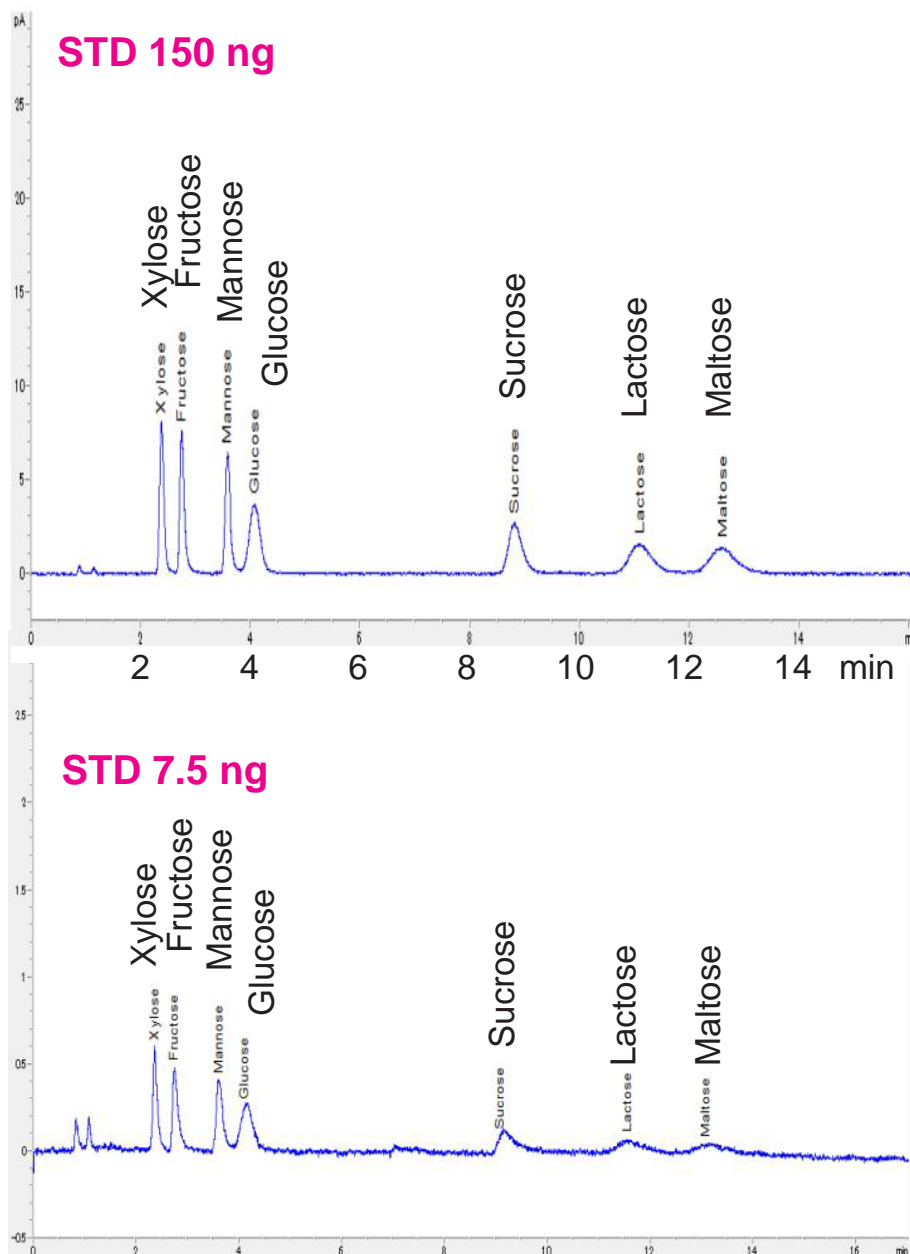

Unison UK-Amino, 250 x 3 mm
 acetonitrile / water = 92 / 8
 1.5 mL/min
 Pre-column temp.: 65 °C, Post-column temp.: 30 °C
 Corona ultra (Filter: High, Nebulizer Temp.: OFF)
 1.5 µL (150ng or 7.5ng)

Courtesy of Kyoko Fukushima, LMS CO., LTD. / JAPAN