


Cadenza CX-C18

50 x 2 mm

Application

LC-MS: Food preservatives

食品保存料のLC-MS分析

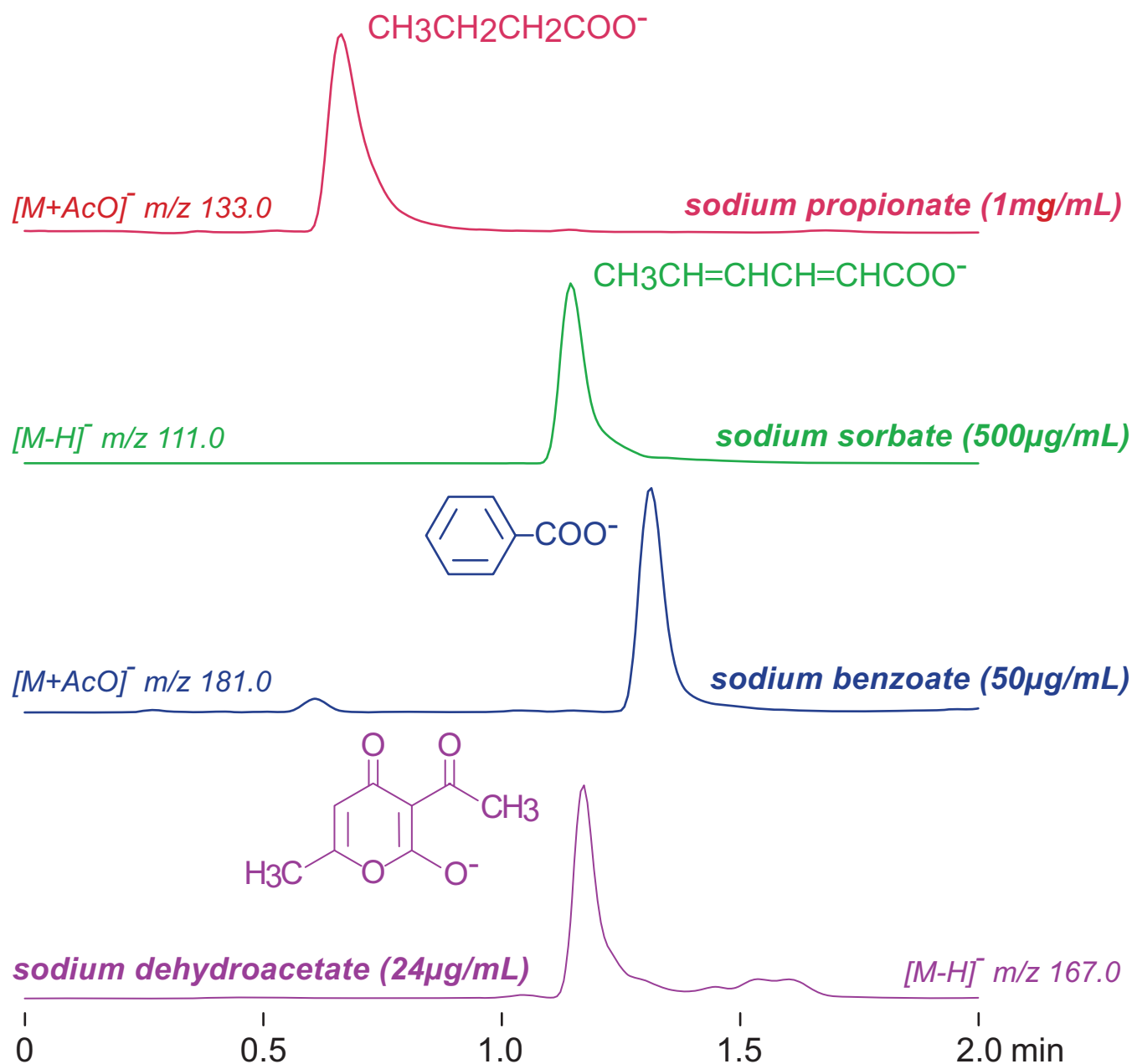

Cadenza CX-C18, 50 x 2 mm  
 A: water / acetic acid = 100 / 0.1  
 B: acetonitrile / acetic acid = 100 / 0.1  
 20-90%B (0-1.2min), 90%B (1.2-2min)  
 0.3mL/min (7MPa), 37°C, MS (ESI, SIM, Negative)  
 5 $\mu\text{L}$  (0.1% AcOH in 15% ACN)