




Unison UK-Phenyl

150 x 3 mm

Application

LC-MS: Catechins in Black Tea with tea bag

ティーバッグ紅茶中のカテキン類のLC-MS分析

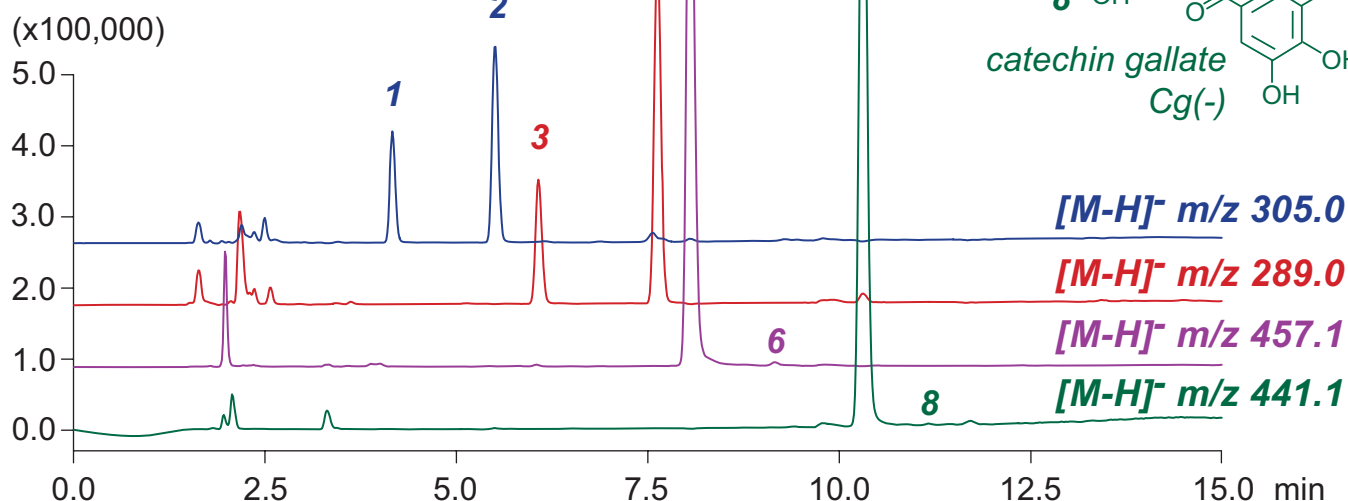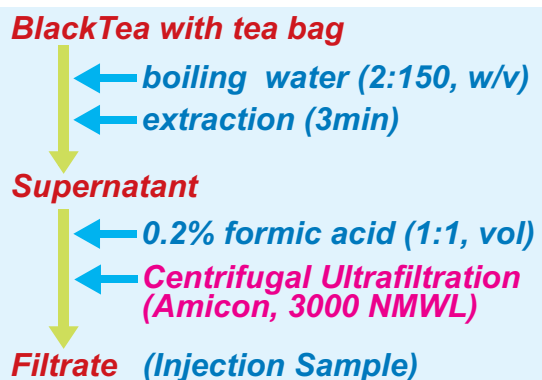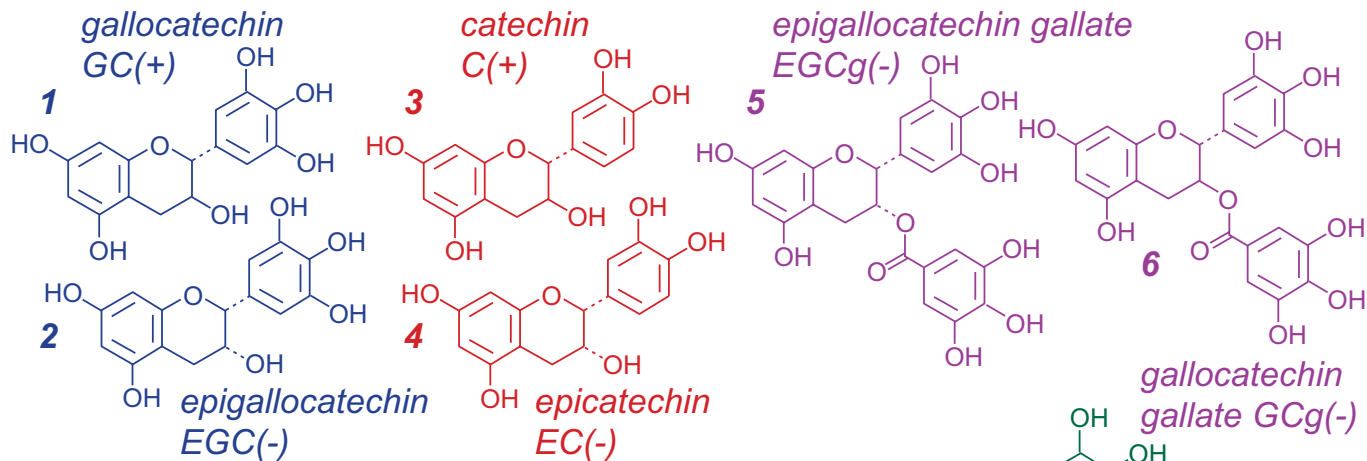

Unison UK-Phenyl, 150 x 3 mm
 A: water / formic acid = 100 / 0.1
 B: methanol / water / formic acid = 40 / 60 / 0.1
 20-100%B (0-12min), 100%B (12-15min)
 0.4mL/min (8MPa), 37°C, 5µL
 MS (ESI, SIM, Negative)