


Scherzo SW-C18  
 Scherzo SM-C18  
 Scherzo SS-C18

100 x 4.6 mm

HPLC COLUMN APPLICATION

## Mesalazine (Anti-inflammatory) and isomers

メサラジン（抗炎症薬）と異性体

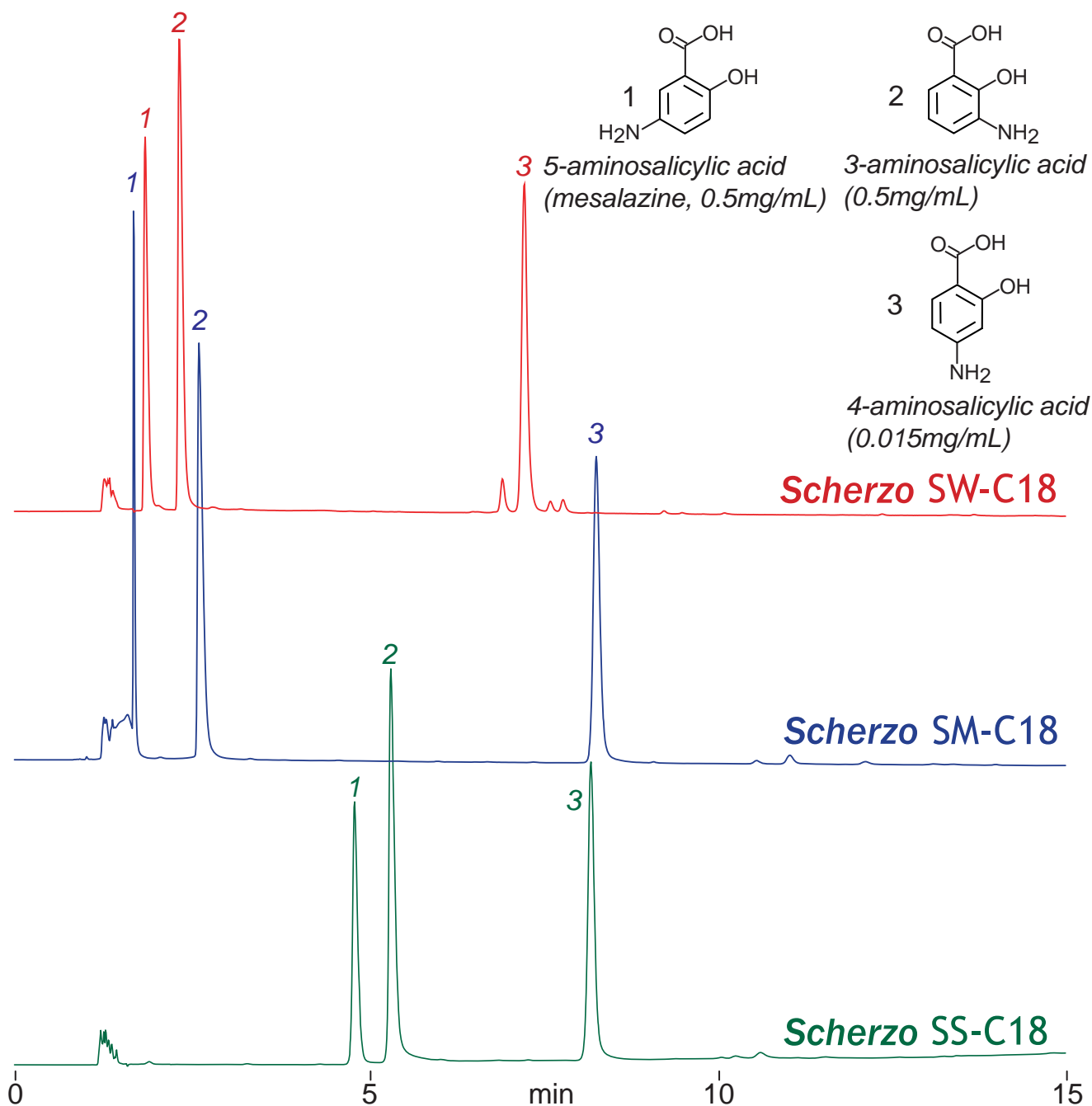

100 x 4.6 mm

A: water / formic acid = 100 / 0.2

B: 100mM ammonium acetate / acetonitrile = 70 / 30

10-75%B (0-15min)

1.0 mL/min (12MPa), 37°C, 260nm

10µL (0.015-0.5mg/mL, water / DMSO / HCOOH = 50 / 50 / 1)