




Scherzo SW-C18

100 x 3 mm

HPLC COLUMN APPLICATION

LC-MS: Styrene urinary metabolites

尿中スチレン代謝物の LC-MS 分析

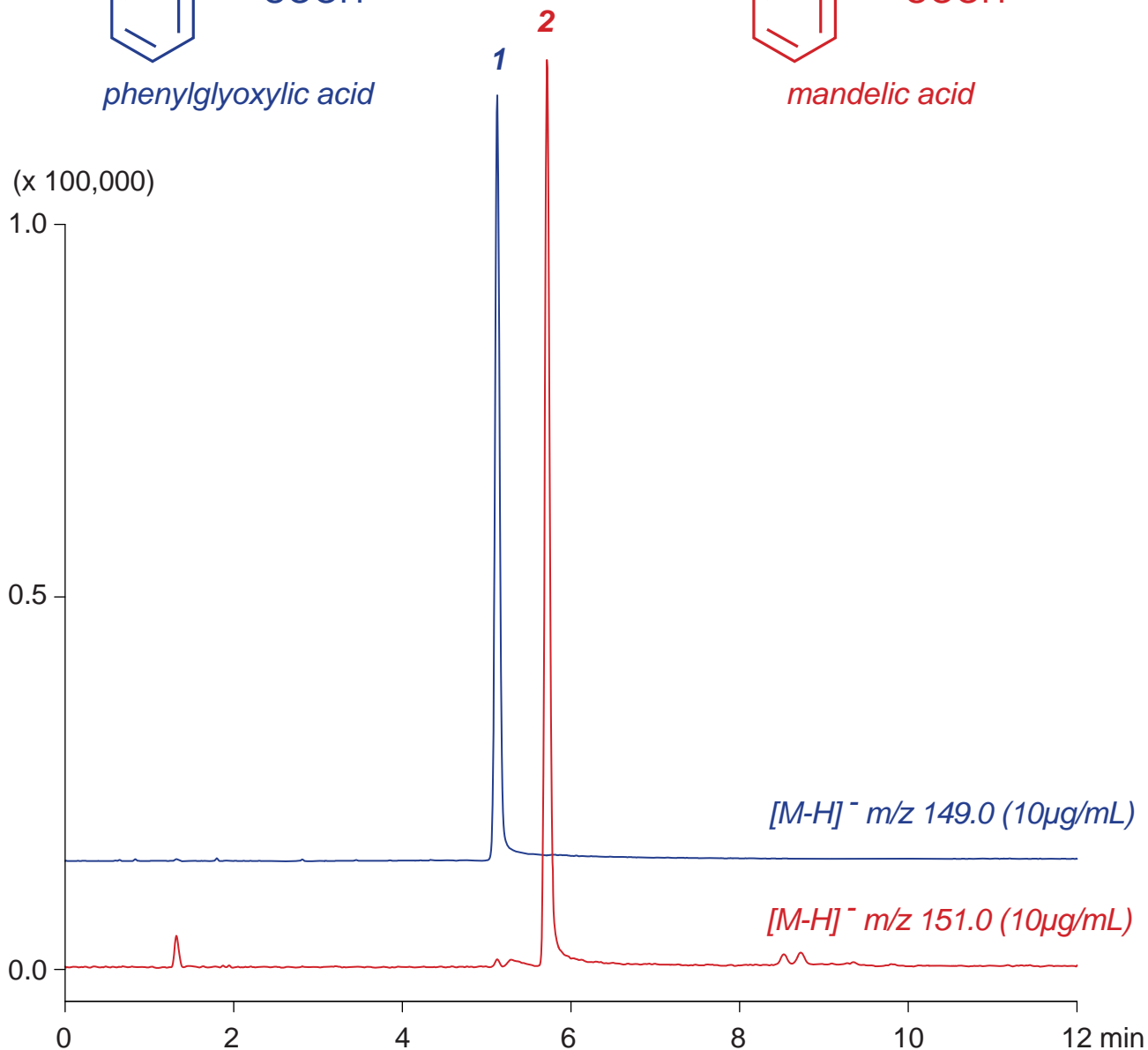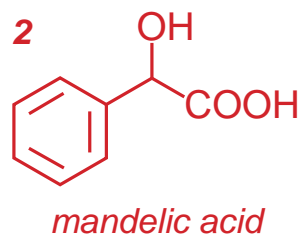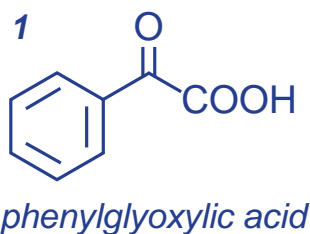

Scherzo SW-C18, 100 x 3 mm

A: water / formic acid = 100 / 0.5

B: acetonitrile / 100mM ammonium formate = 50 / 50

0-100%B (0-10min), 100%B (10-12min)

0.4mL/min (7MPa), 37°C, Single Quad. MS (ESI)

5µL (10µg/mL, 1% formic acid)