



Unison UK-Phenyl  
Unison UK-C8  
Unison UK-C18

Cadenza CX-C18  
Cadenza CW-C18  
Cadenza CD-C18 100 x 4.6 mm

HPLC COLUMN APPLICATION

## Rhododendrol (Tyrosinase inhibitor)

ロドデンドロール (チロシナーゼ阻害剤)

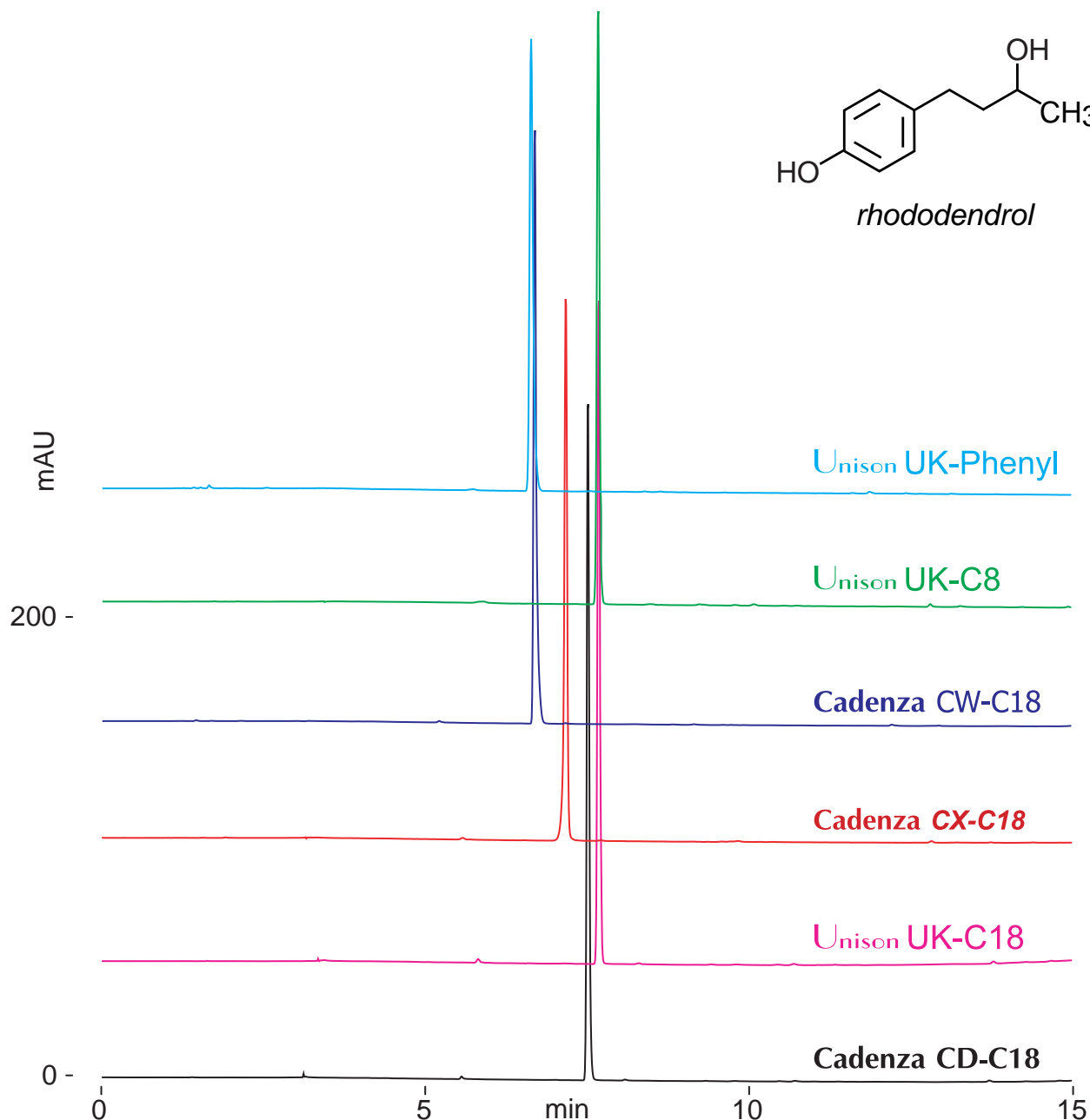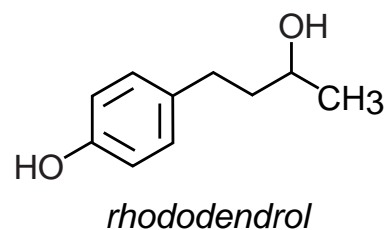

100 x 4.6 mm

A: 100mM ammonium formate

B: 100mM ammonium formate / acetonitrile = 10 / 90

0-90%B (0-15min)

1.0mL/min (9MPa), 37°C, 280nm

10µL (0.3mg/mL, 100mM ammonium formate)