

Unison UK-Phenyl Cadenza CX-C18



Unison UK-C8 Cadenza CW-C18

Unison UK-C18 Cadenza CD-C18 100 x 4.6 mm

HPLC COLUMN APPLICATION

Piceid (Resveratrol derivatives)

ピセイド (レスベラトロール誘導体)

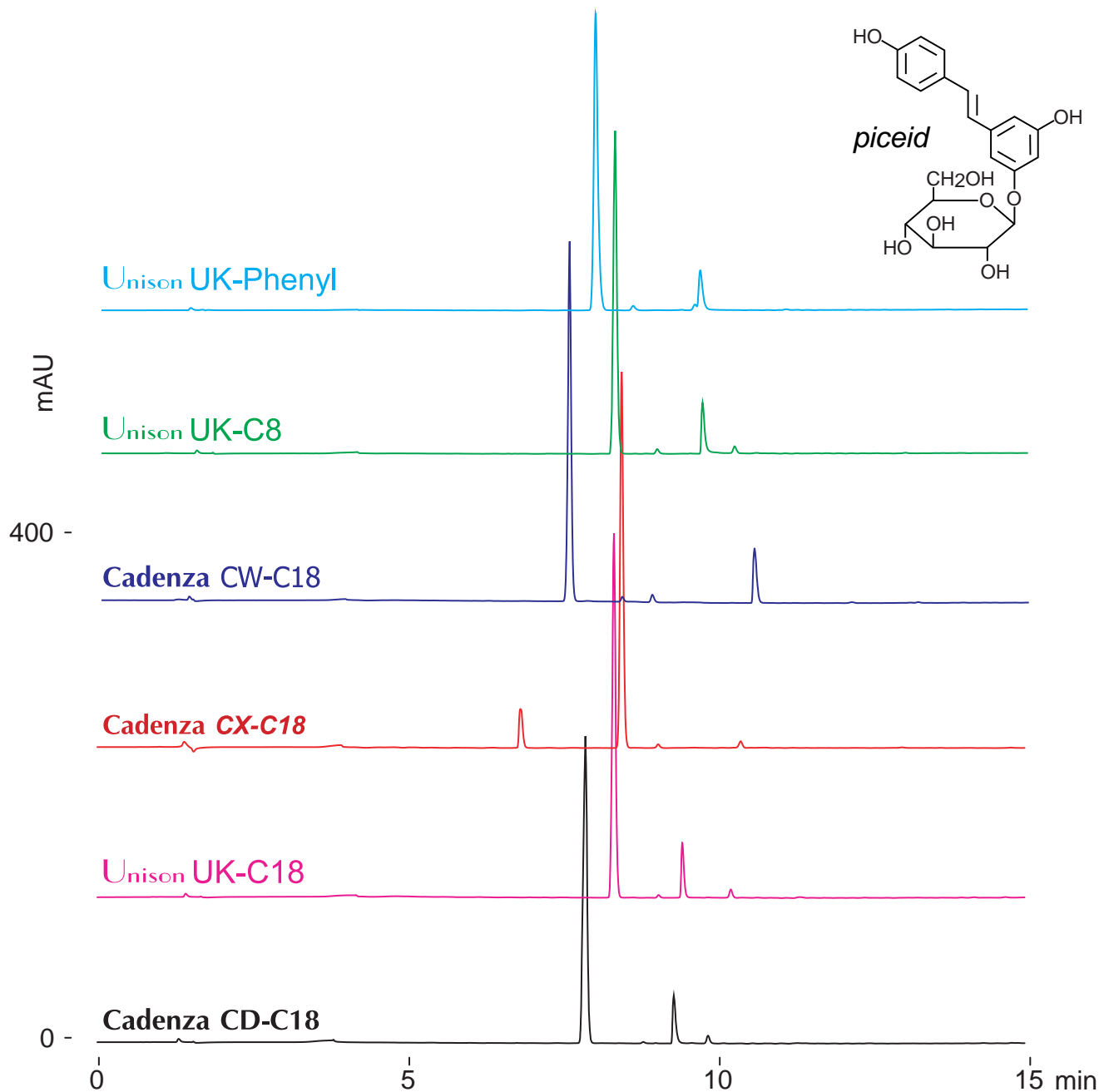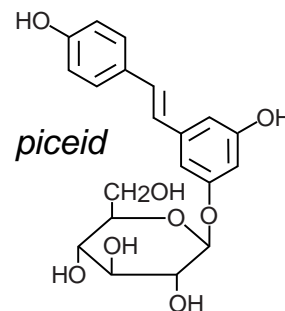

100 x 4.6 mm

A: water / formic acid = 100 / 0.1

B: water / acetonitrile / formic acid = 10 / 90 / 0.1

0%B (0-1min), 0-70%B (1-15min)

1.0mL/min (9MPa), 37°C, 260nm

10µL (0.4mg/mL, water / acetonitrile / formic acid = 50 / 50 / 1)