

Unison UK-Phenyl Cadenza CX-C18



Unison UK-C8 Cadenza CW-C18

Unison UK-C18 Cadenza CD-C18 100 x 4.6 mm

HPLC COLUMN APPLICATION

Aloe emodin (Stimulant laxative)

アロエエモジン (刺激性下剤)

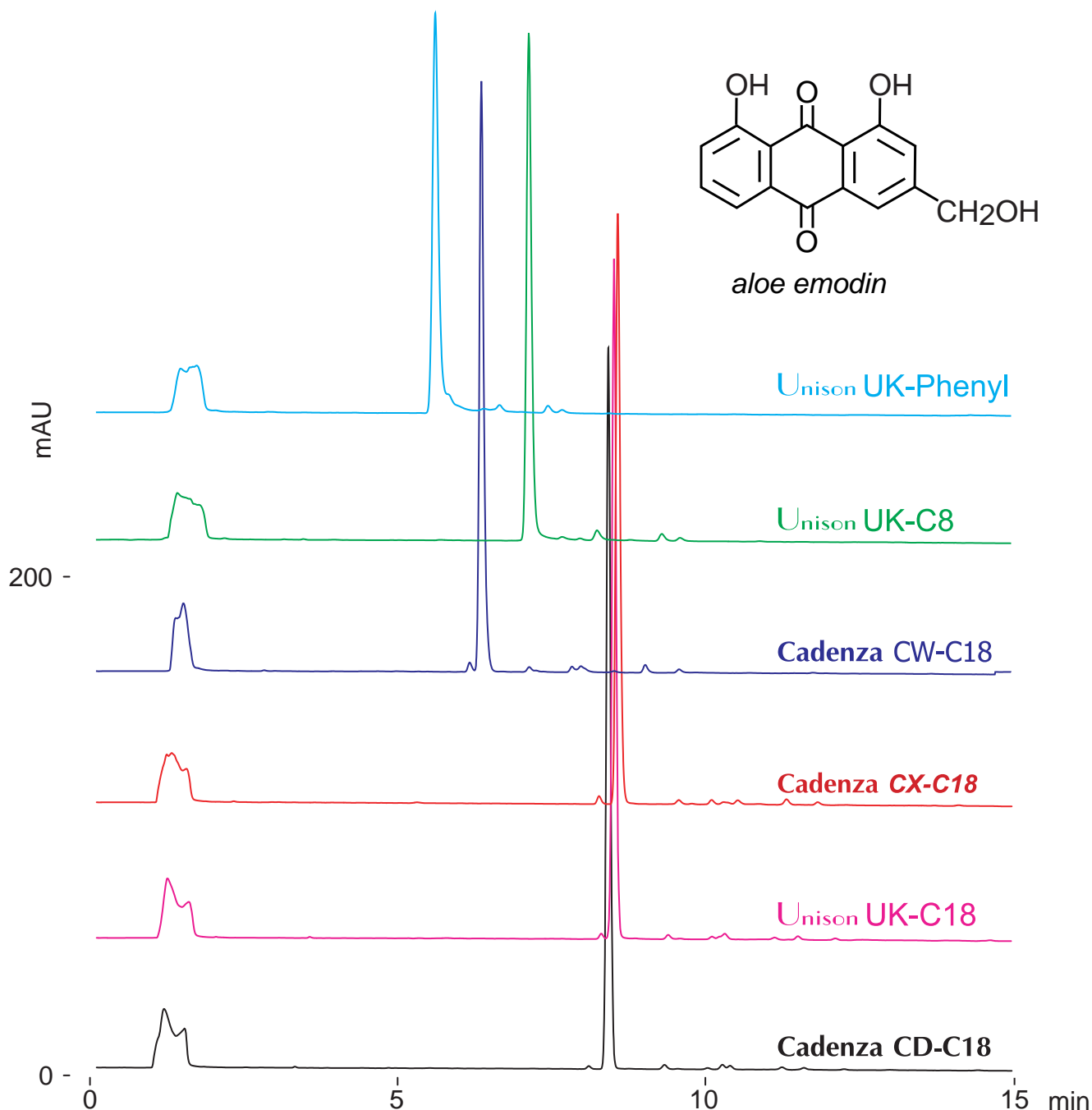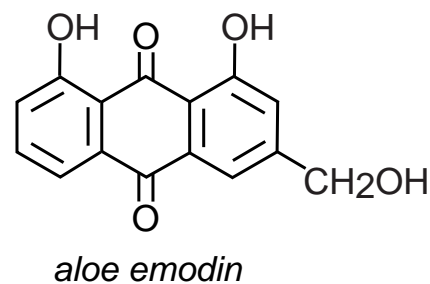

100 x 4.6 mm

A: 100mM ammonium formate / acetonitrile = 90 / 10

B: 100mM ammonium formate / acetonitrile = 10 / 90

20-100%B (0-15min)

1.0mL/min (10MPa), 37°C, 260nm

10µL (0.075mg/mL, 100mM ammonium formate / DMF = 10 / 90)