




Cadenza CD-C18

100 x 3 mm

HPLC COLUMN APPLICATION

LC-MS: Capsaicin (Capsaicinoid)

カプサイシン (カプサイシノイド) の LC-MS 分析

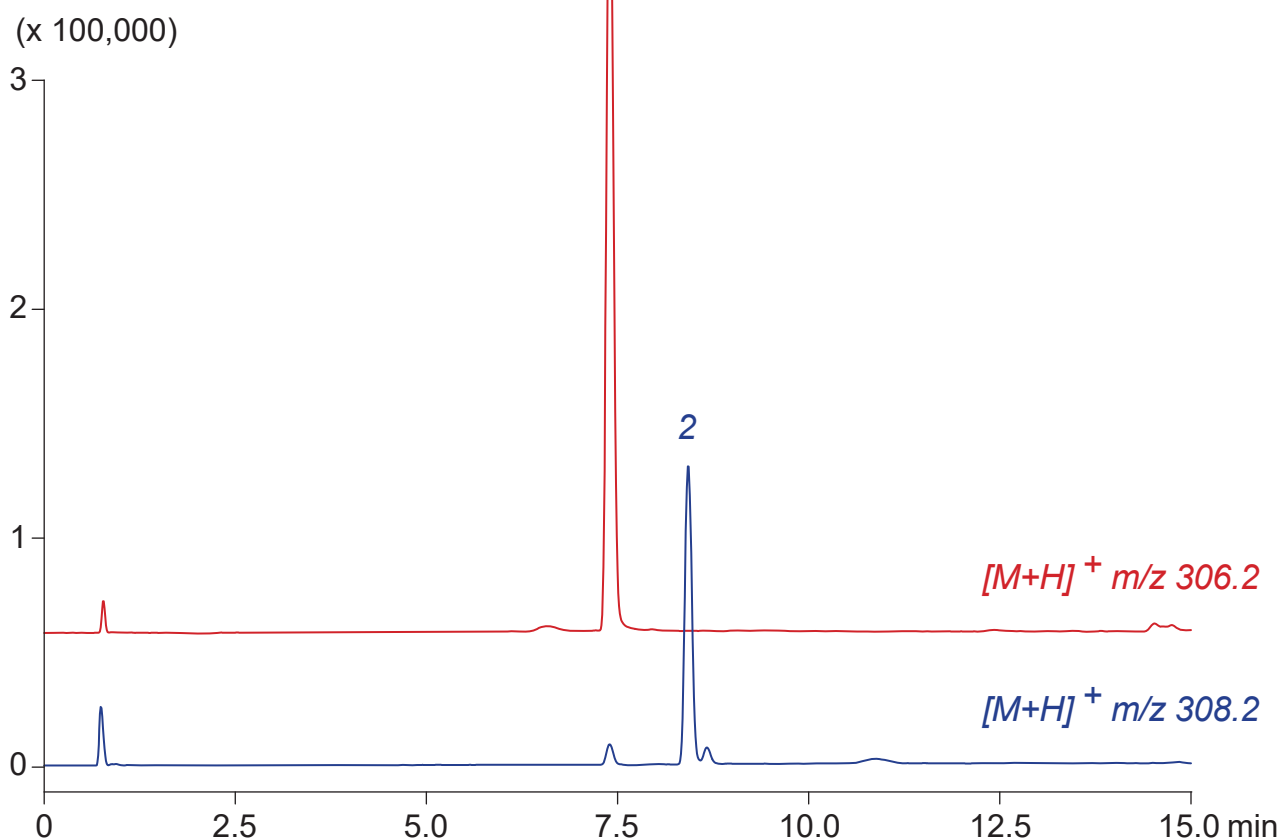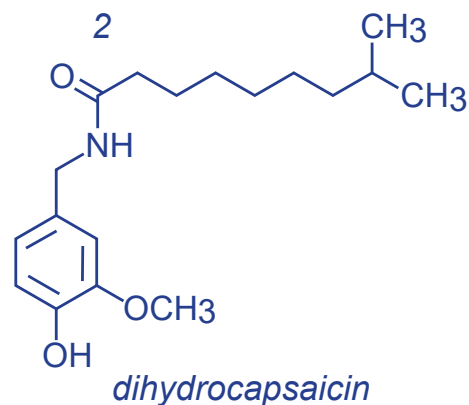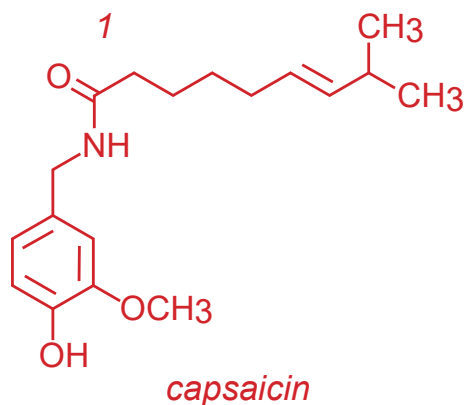

Cadenza CD-C18, 100 x 3 mm

A: acetonitrile / water / formic acid = 10 / 90 / 0.1

B: acetonitrile / water / formic acid = 90 / 10 / 0.1

30-100%B (0-12min), 100%B (12-15min)

0.4mL/min (7MPa), 37°C, Single Quad. MS (ESI)

5µL (1µg/mL, ACN / water / HCOOH = 50 / 50 / 1)