




**Scherzo SM-C18 MF**

100 x 2 mm

**HPLC COLUMN APPLICATION**

**LC-MS/MS: Glyphosate, Glufosinate**

グリホサート, グルホシネートの LC-MS/MS 分析

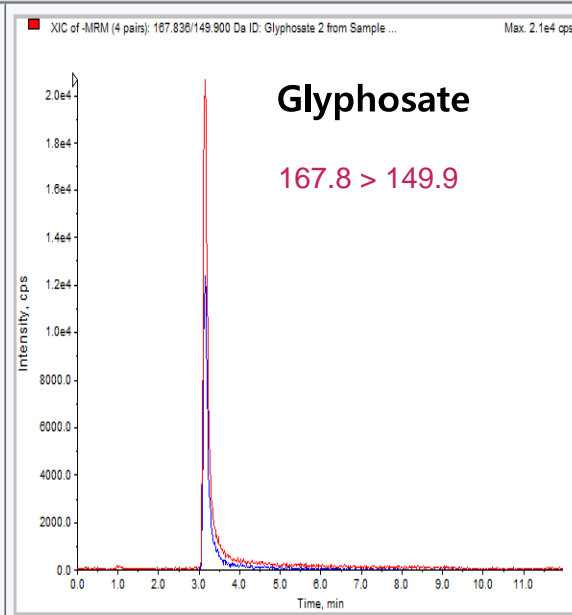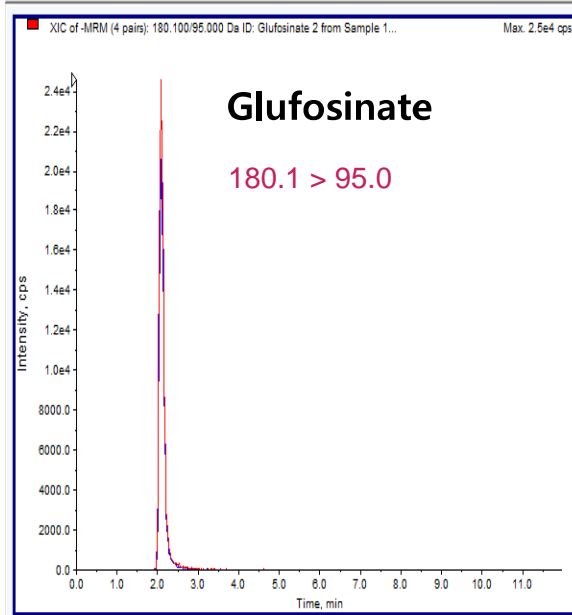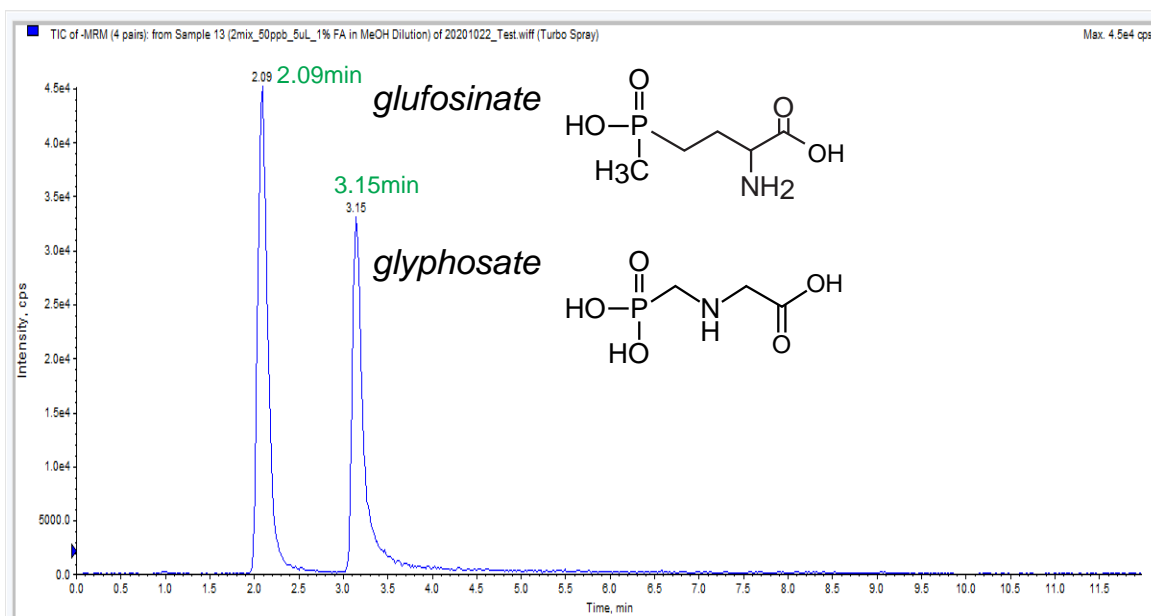

Scherzo SM-C18 MF, 100 x 2 mm

A: 0.1% formic acid in 20% methanol

B: 1% formic acid in 20% methanol

0%B (0-1min), 0-100%B (1-4min), 100%B (4-7min),

100-0%B (7-7.1min), 0%B (7.1-15min)

0.2mL/min, 40°C, Triple Quad. MS (ESI, Negative)

5µL (50ppb)

Courtesy of PEAKMAN SP, KOREA