


Cadenza CD-C18

50 x 2 mm

HPLC COLUMN APPLICATION

LC-MS/MS: Spinosad

スピノサドの LC-MS/MS 分析

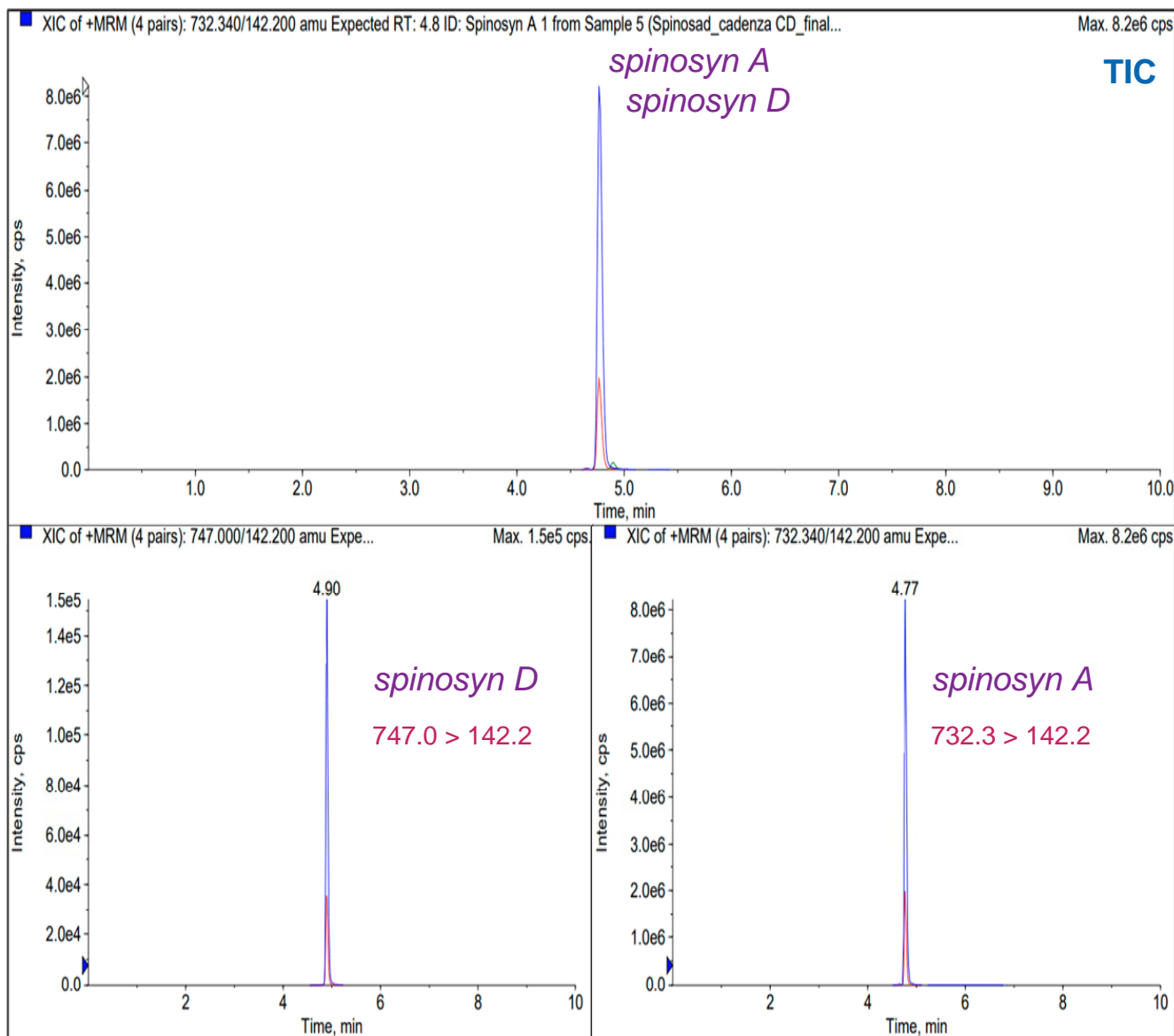

Cadenza CD-C18, 50 x 2 mm

A: 0.1% formic acid in water

B: 0.1% formic acid + 5mM ammonium acetate in methanol

10-100%B (0-4min), 100%B (4-5min), 100-10%B (5-7min), 10%B (7-10min)

0.3mL/min, 40°C, Triple Quad. MS (ESI, Positive)

5µL

Courtesy of PEAKMAN SP, KOREA