




Unison UK-C18

100 x 2 mm

HPLC COLUMN APPLICATION

LC-MS/MS: Stevia sweeteners

ステビア甘味料の LC-MS/MS 分析

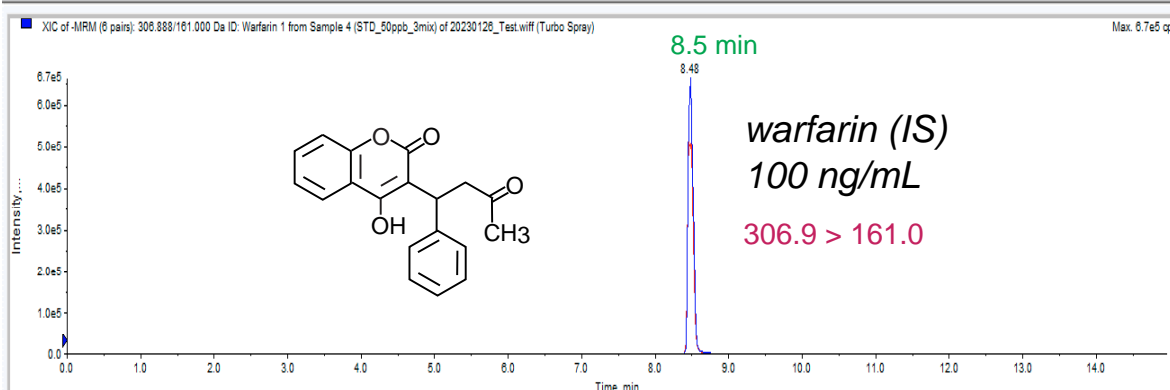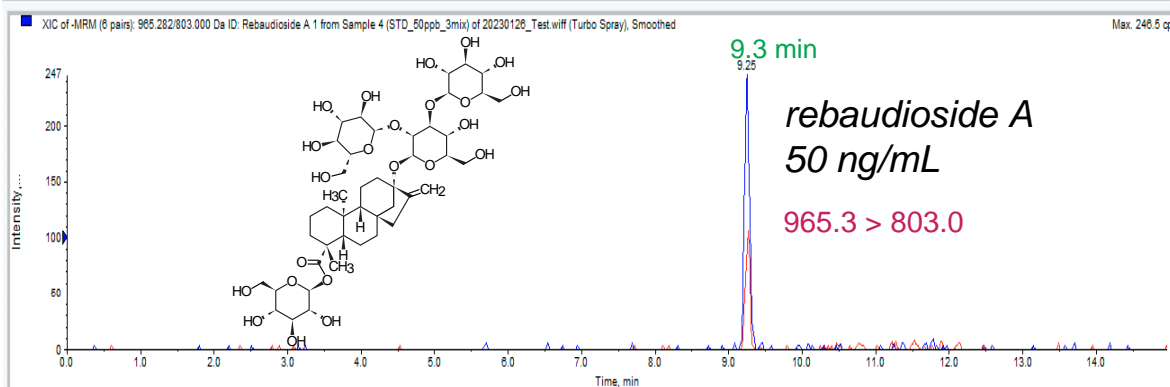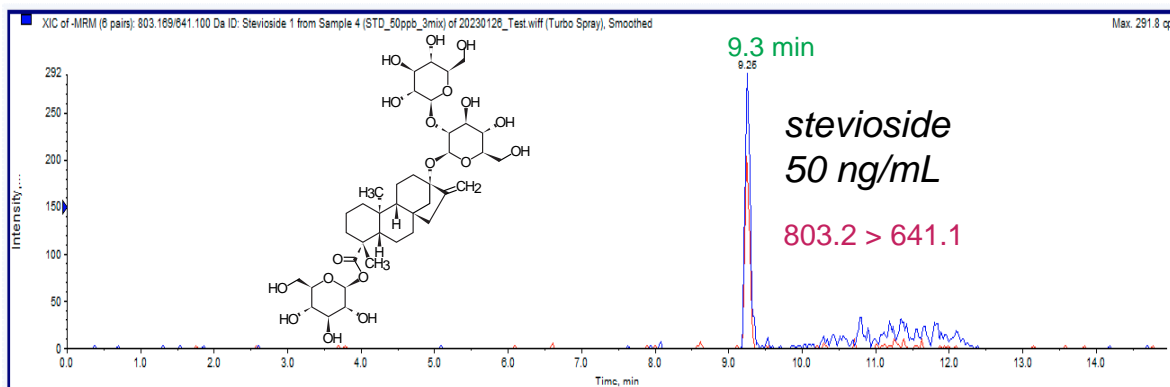

Unison UK-C18, 100 x 2 mm

A: 10mM ammonium acetate in water

B: 10mM ammonium acetate in methanol

5%B (0-3min), 5-95%B (3-7min), 95%B (7-10min),

95-5%B (10-10.1min), 5%B (10.1-15min)

0.2mL/min, 40°C, Triple Quad. MS (ESI, Negative)

5µL

Courtesy of PEAKMAN SP, KOREA