




Scherzo SS-C18

150 x 3 mm

HPLC COLUMN APPLICATION

LC-MS/MS: Chlorate

塩素酸の LC-MS/MS 分析

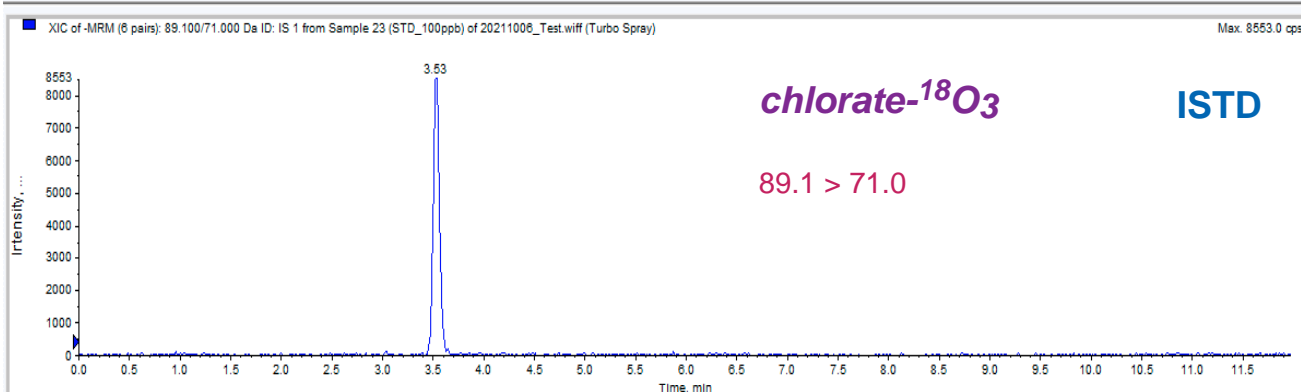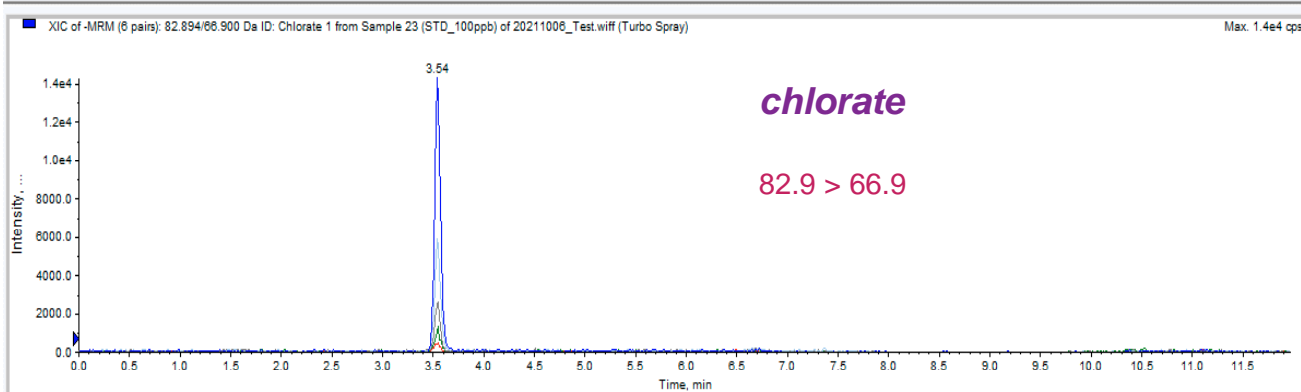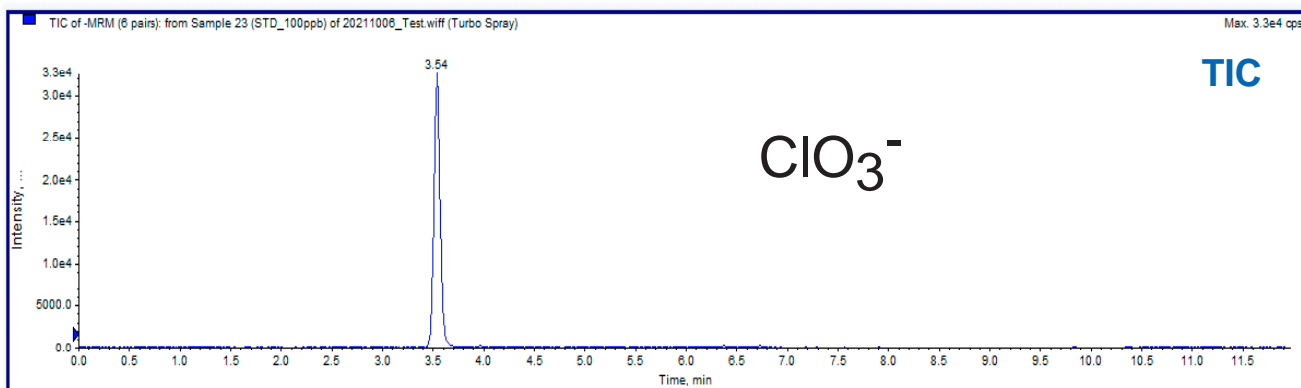

Scherzo SS-C18, 150 x 3 mm

A: 50mM ammonium formate + 0.45% formic acid in 40% acetonitrile

B: acetonitrile

0%B (0-4min), 0-90%B (4-6min), 90%B (6-8min),

90-0%B (8-8.1min), 0%B (8.1-12min)

0.4mL/min, 35°C, Triple Quad. MS (ESI, Negative)

5µL (100 ppb)

Courtesy of PEAKMAN SP, KOREA