


Unison UK-C18 HT 100 x 2 mm HPLC COLUMN APPLICATION

LC-MS/MS: Antibiotics (7)
抗生物質の LC-MS/MS 分析 (7)

Antibiotics 107mix 10min method

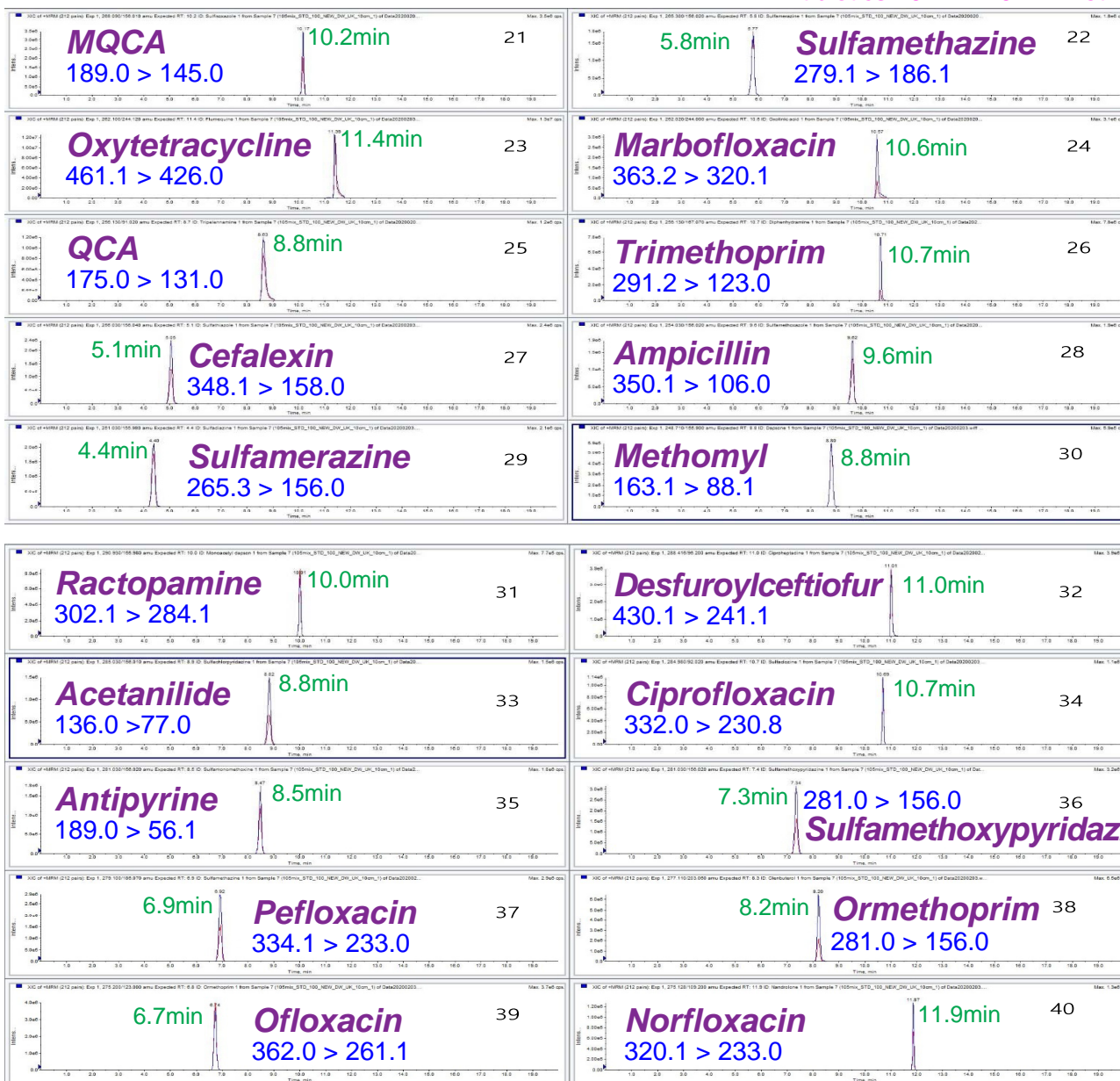

Unison UK-C18 HT, 100 x 2 mm
 A: 0.1% formic acid in water
 B: 0.1% formic acid in acetonitrile
 10%B (0-0.5min), 10-15%B (0.5-3min), 15-95%B (3-6min),
 95%B (6-8min), 95-10%B (8-8.1min), 10%B (8.1-10min)
 0.4mL/min, 40°C, Triple Quad. MS (ESI, Positive/Negative)
 5µL

Courtesy of PEAKMAN SP, KOREA