







Unison UK-C18 HT

100 x 2 mm

HPLC COLUMN APPLICATION

LC-MS/MS: Anticoccidial agents

抗コキシジウム薬の LC-MS/MS 分析

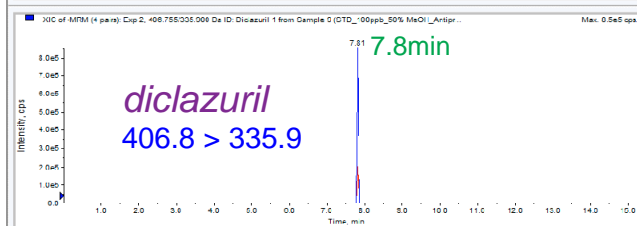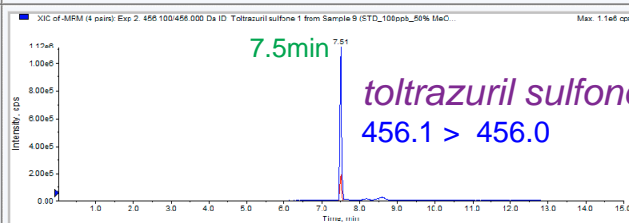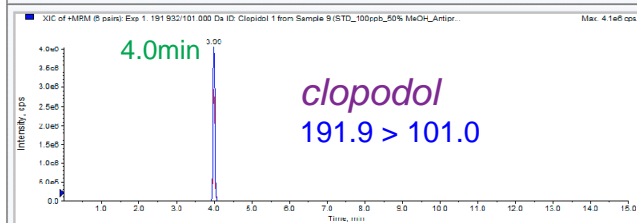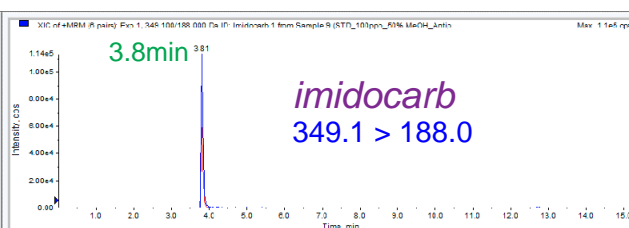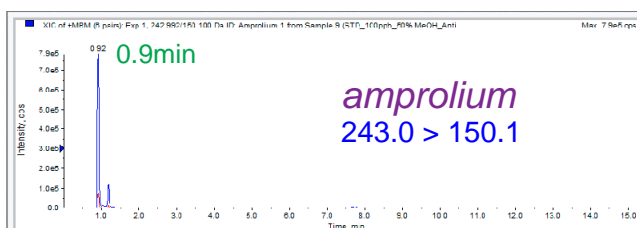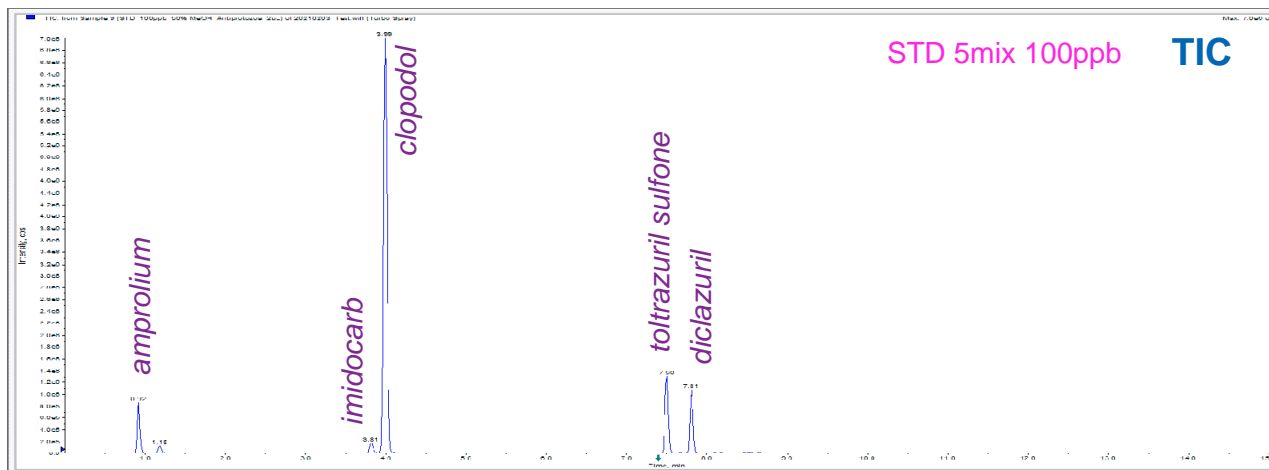

Unison UK-C18 HT, 100 x 2 mm
 A: 0.1% formic acid in water
 B: 0.1% formic acid in acetonitrile
 5%B (0-1min), 5-95%B (1-8min), 95%B (8-11min),
 95-5%B (11-11.1min), 5%B (11.1-15min)
 0.3mL/min, 40°C, Triple Quad. MS (ESI, Positive/Negative)
 2µL (100 ppb)

Courtesy of PEAKMAN SP, KOREA