








Cadenza CD-C18 HT

150 x 2 mm

HPLC COLUMN APPLICATION

LC-MS/MS: Aminoglycoside antibiotics (2)

アミノグリコシド系抗生物質の LC-MS/MS 分析 (2)

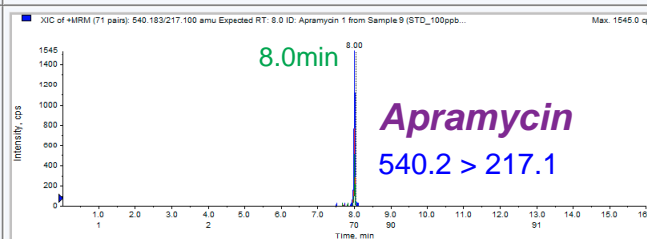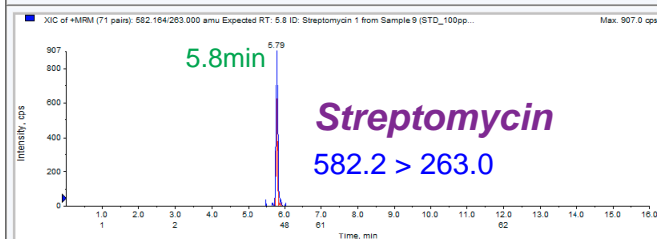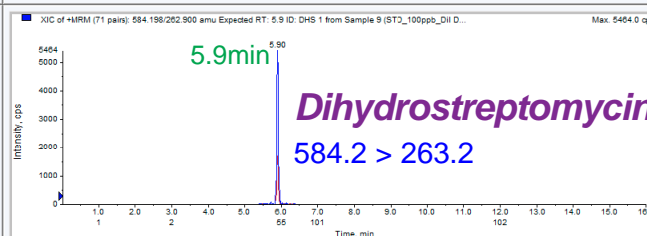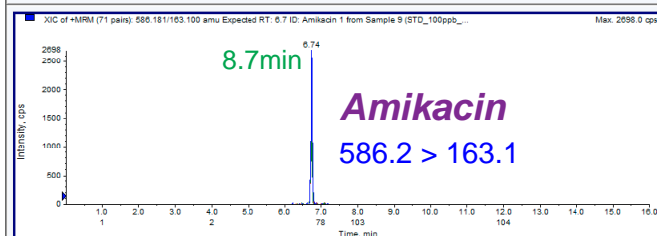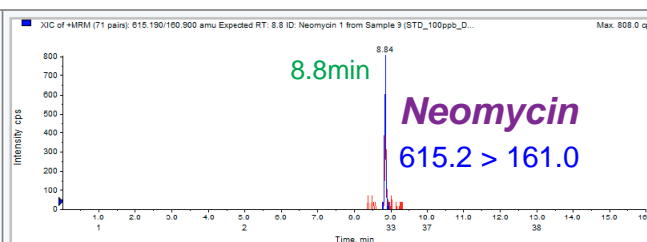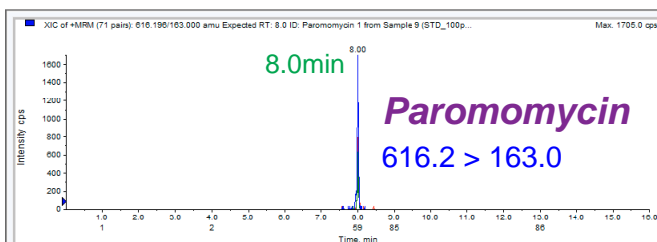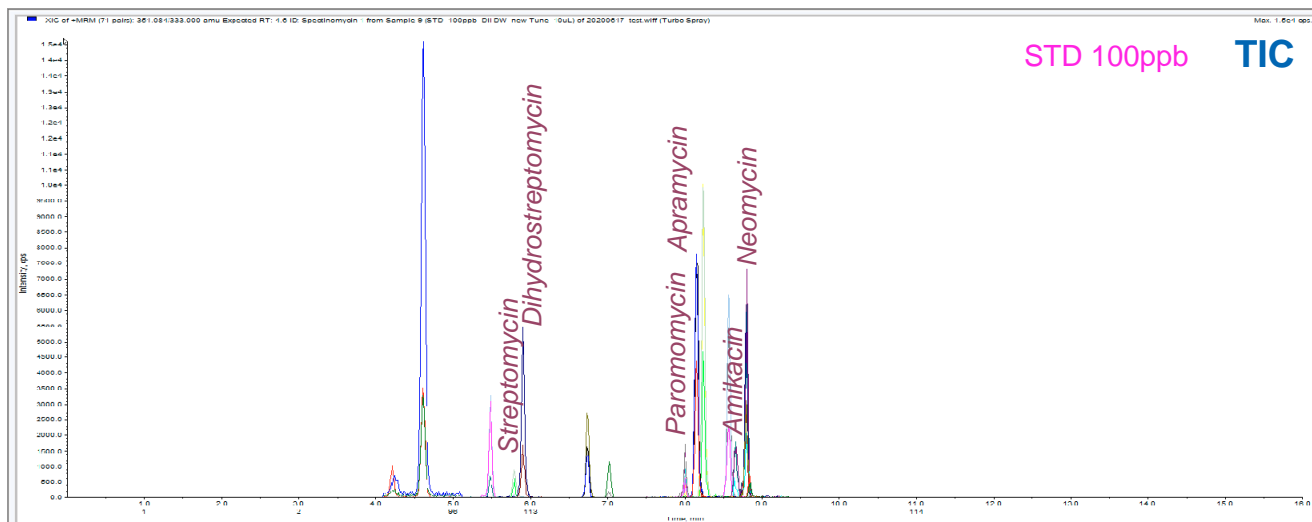

Cadenza CD-C18 HT, 150 x 2 mm
 A: 20mM HFBA in 5% acetonitrile
 B: 20mM HFBA in 50% acetonitrile
 20%B (0-0.5min), 20-95%B (0.5-9min), 95%B (9-12min),
 95-20%B (12-12.1min), 20%B (12.1-16min)
 0.3mL/min, 40°C, Triple Quad. MS (ESI, Positive)
 10µL (100 ppb)

Courtesy of PEAKMAN SP, KOREA