


Unison UK-Phenyl Cadenza CX-C18
 Unison UK-C8 Cadenza CW-C18
 Unison UK-C18 Cadenza CD-C18

100 x 4.6 mm

HPLC COLUMN APPLICATION

Monocyclic terpene isomers

単環式モノテルペン異性体

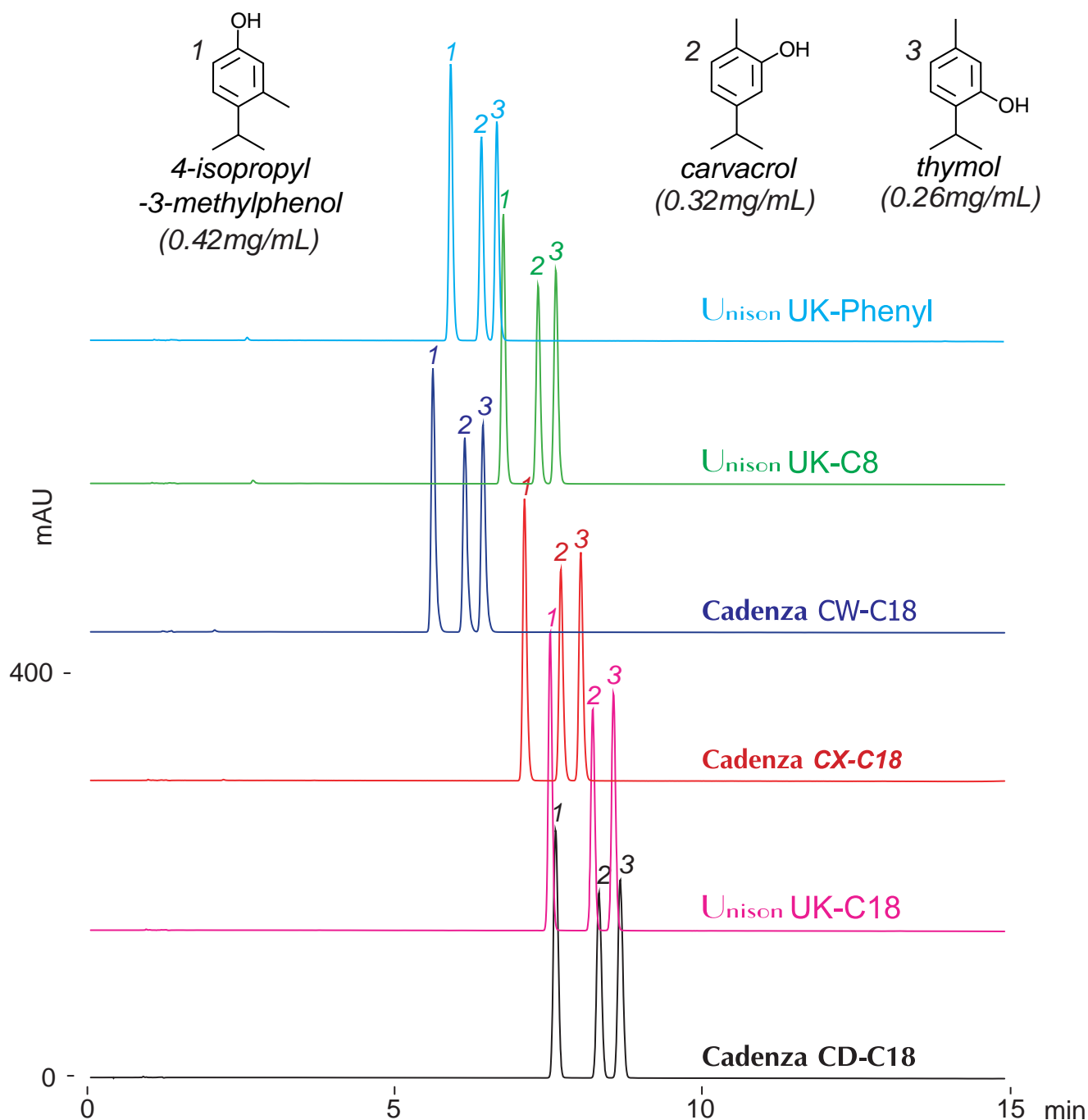

100 x 4.6 mm

A: 100mM ammonium formate / acetonitrile = 90 / 10

B: 100mM ammonium formate / acetonitrile = 10 / 90

30-100%B (0-15min)

1.0mL/min (10MPa), 37°C, 280nm

10µL (0.26-0.42mg/mL, 100mM ammonium formate / acetonitrile = 80 / 20)