

Unison UK-Phenyl

Unison UK-C8

Unison UK-C18

Cadenza CW-C18


Cadenza CD-C18

100 x 4.6 mm


HPLC COLUMN APPLICATION

Organic Redox-Flow Battery positive electrolyte ((Ferrocenylmethyl)trimethylammonium Chloride)

有機レドックスフロー電池正極電解質
((フェロセニルメチル)トリメチルアンモニウムクロリド)

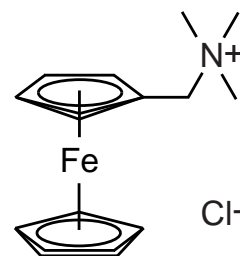

(ferrocenylmethyl)trimethylammonium chloride

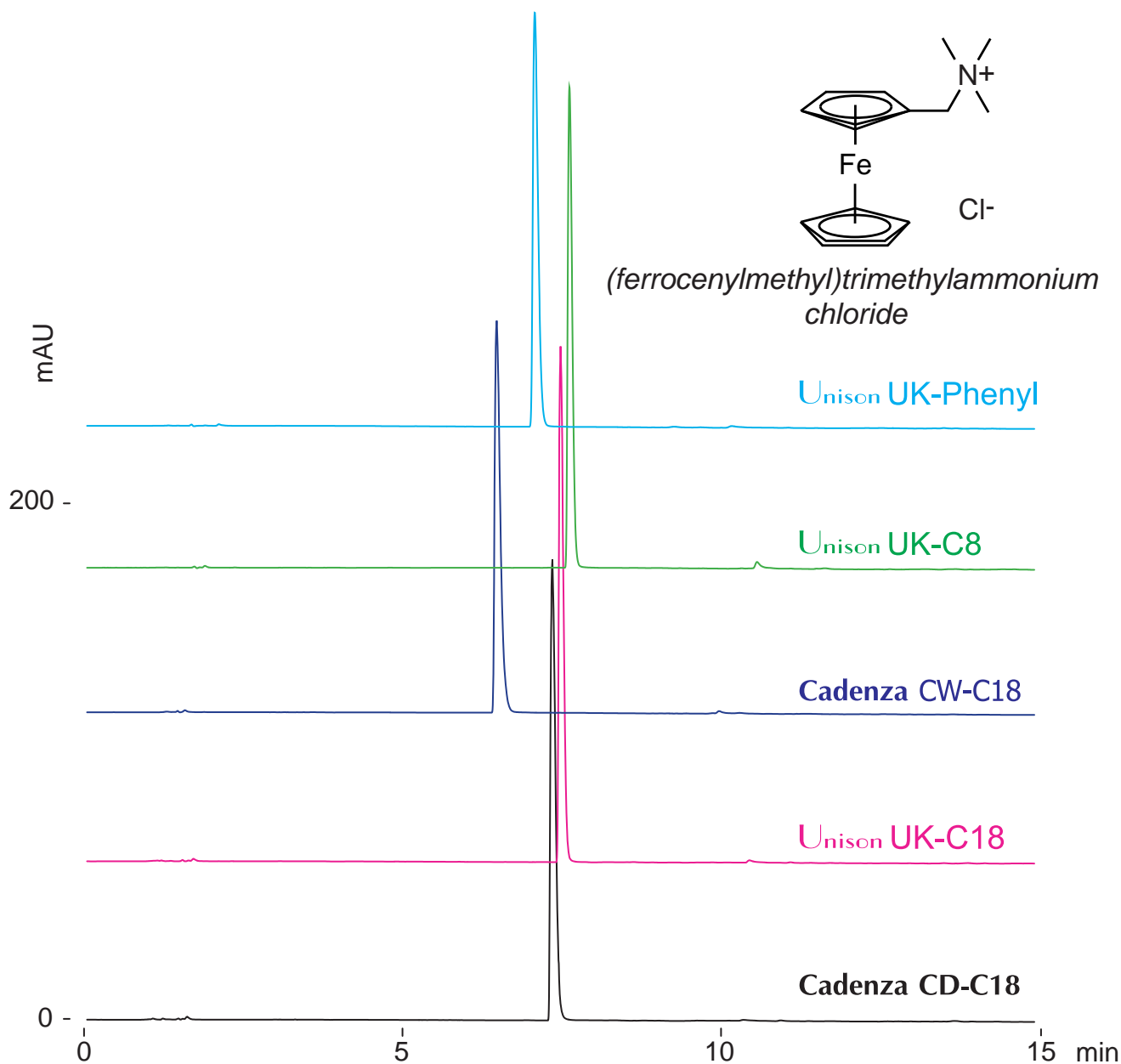

100 x 4.6 mm

A: 100mM ammonium formate / acetonitrile = 90 / 10

B: 100mM ammonium formate / acetonitrile = 10 / 90

0%B (0-1min), 0-60%B (1-15min)

1.0mL/min (9MPa), 37°C, 260nm

10µL (0.25mg/mL, 100mM ammonium formate)