


Cadenza CD-C18

100 x 4.6 mm


HPLC COLUMN APPLICATION

Organic Redox-Flow Battery positive electrolyte  
(TEMPO Free Radical)


有機レドックスフロー電池正極電解質 (TEMPOフリーラジカル)

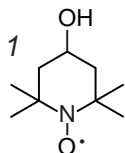


4-hydroxy-2,2,6,6-tetramethyl piperidine 1-oxyl free radical

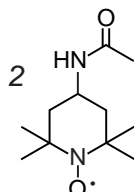


4-acetamido-2,2,6,6-tetramethyl piperidine 1-oxyl free radical

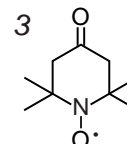


4-oxo-2,2,6,6-tetramethyl piperidine 1-oxyl free radical

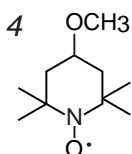

4-methoxy-2,2,6,6-tetramethyl piperidine 1-oxyl free radical

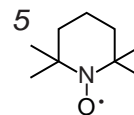

2,2,6,6-tetramethyl piperidine 1-oxyl free radical

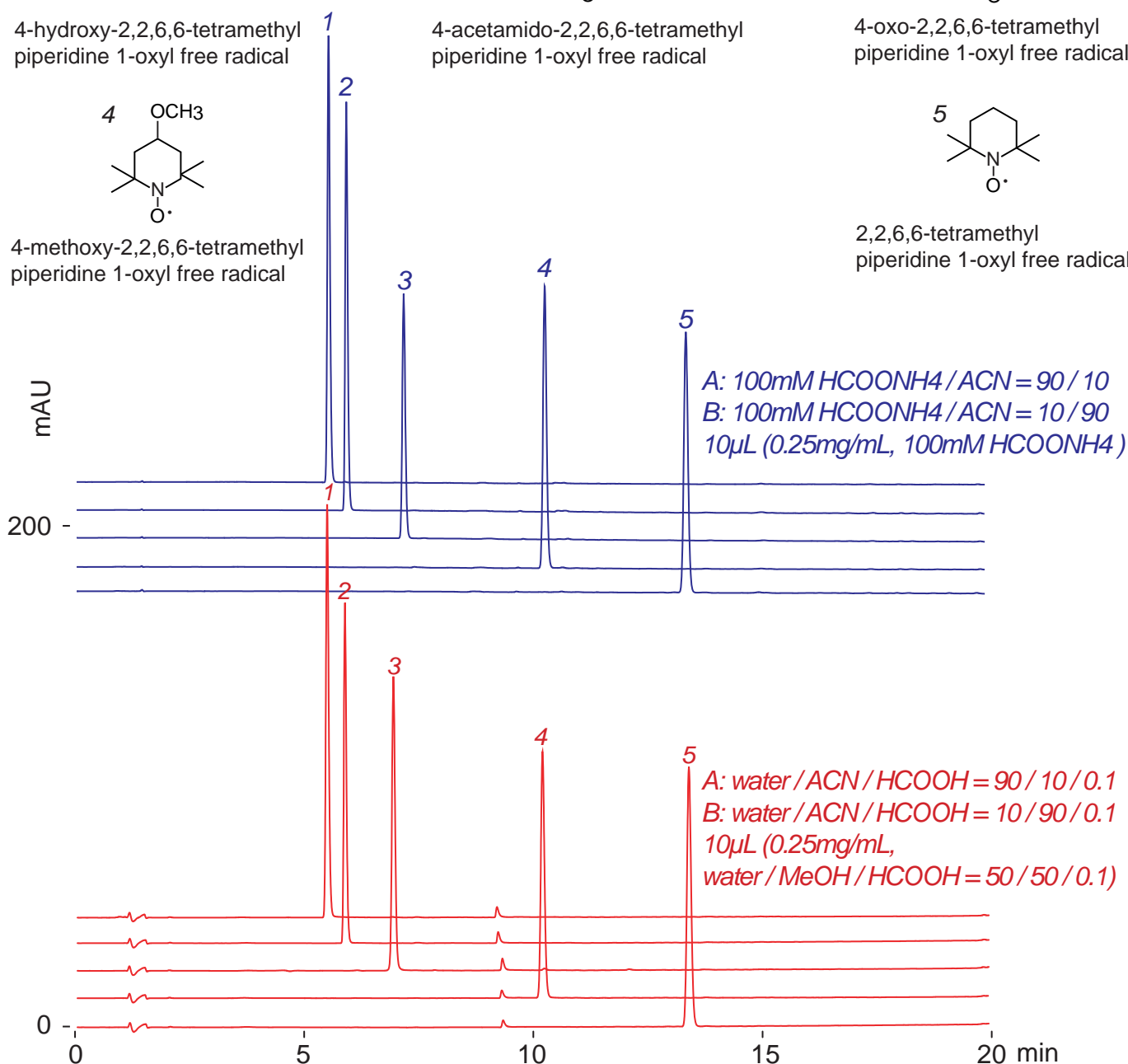

Cadenza CD-C18, 100 x 4.6 mm  
0%B (0-1min), 0-80%B (1-20min)  
1.0mL/min (9MPa), 37°C, 260nm